
autoCorrelator

Release 1.0

CAS

Jun 01, 2021

1	Introduction	3
2	Calibration	7
3	Device Scenes	9
4	Troubleshooting	11
5	Indices and tables	13

Contents:

Introduction

The AutoCorrelator device is designed to provide an online determination of the pulse duration using a single-shot auto correlator ¹.

The measurement of the time profile of pulses is based on the following principle, graphically displayed in Fig. 1.1. The input beam is sent to a beam-splitter; the two identical beams propagate along two distinct optical paths until they intersect in a non-linear crystal. Here, due to the high-intensity of the beams, a second harmonic beam (SH) is created and its integrated energy is measured by a CCD camera located after the crystal.

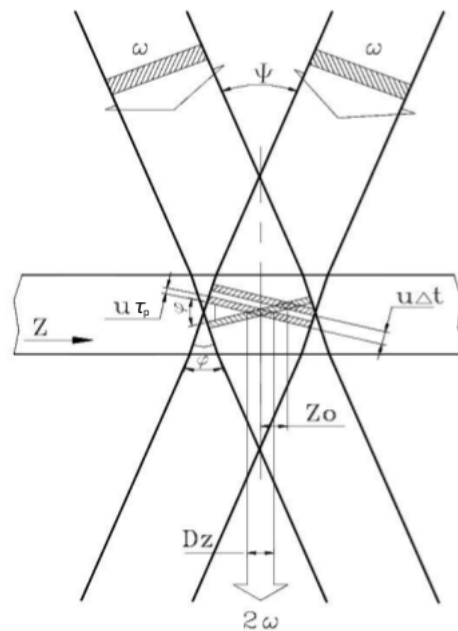


Fig. 1.1: The diagram describes geometrically the intersection of two identical beams in a crystal and the generation of the second harmonic beam.

The pulse duration of laser pulses can be determined upon measuring the transverse distribution of the energy deposited in the CCD camera. From geometrical considerations in Fig. 1.1, assuming for the incoming beams a rectangular time profile τ_p and uniform transverse intensity profile, it is found that the transverse profile D_z of the second harmonic

¹ RP Photonics Encyclopedia, <https://www.rp-photonics.com/autocorrelators.html>

depends on the pulse duration τ_p of the fundamental beams,

$$D_z = \frac{\tau_p \cdot u}{\sin(\phi/2)}$$

$$\tau_p = D_z \cdot \frac{1}{2} \cdot \frac{\Delta t}{\Delta Z_0}$$

where $u = c/n$ and ϕ are the speed of light and the intersection angle of input beams, respectively, in the crystal with refractive index n . The transverse profile D_z is determined from the data accumulated with the CCD camera available in the system. An example of camera image is presented in Fig. 1.2:

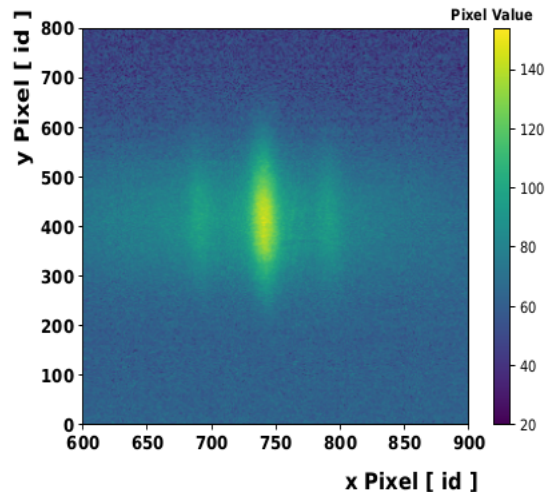


Fig. 1.2: The fundamental beams and the second harmonic beam are detected in the CCD camera located after the non-linear crystal.

The figure shows clearly the deposited energy from the signal of the generated second harmonic beam (central and more intense peak) and of the two fundamental beams (low intensity side signals). The transverse profile D_z is determined as FWHM from the fit to the SH peak.

The angle ϕ cannot be measured with sufficient precision for a reliable extraction of pulse duration τ_p . The way used in the device to determine the pulse duration from the measured transverse profile is presented the calibration section.

The device configuration editor is presented in Fig. 1.3,

The camera device providing the image of the beam profile should be set in the key **input.connectedOutputChannels** of the autocorrelator device. For each camera image the projection along the x-axis is calculated, a fit is performed according to a selectable model (**Beam Shape**) for the time-profile of the pulse, and the peak position and FWHM are determined from the fitting function (**Input Image Peak x-Pos** and **Input Image Peak x-Pos**). The **Fit Error** parameter is an integer flag describing the fit status. If it is equal to 1, 2, 3 or 4, the solution was found, otherwise the solution was not found². The possible fit status values are:

- 0: Improper input parameters were entered,
- 1: The solution converged,
- 2: The number of calls to function has reached default max number,
- 3: Max for relative error is too small, no further improvement in the approximate solution is possible,

² Scipy.org, <https://github.com/scipy/scipy/blob/master/scipy/optimize/minpack.py>

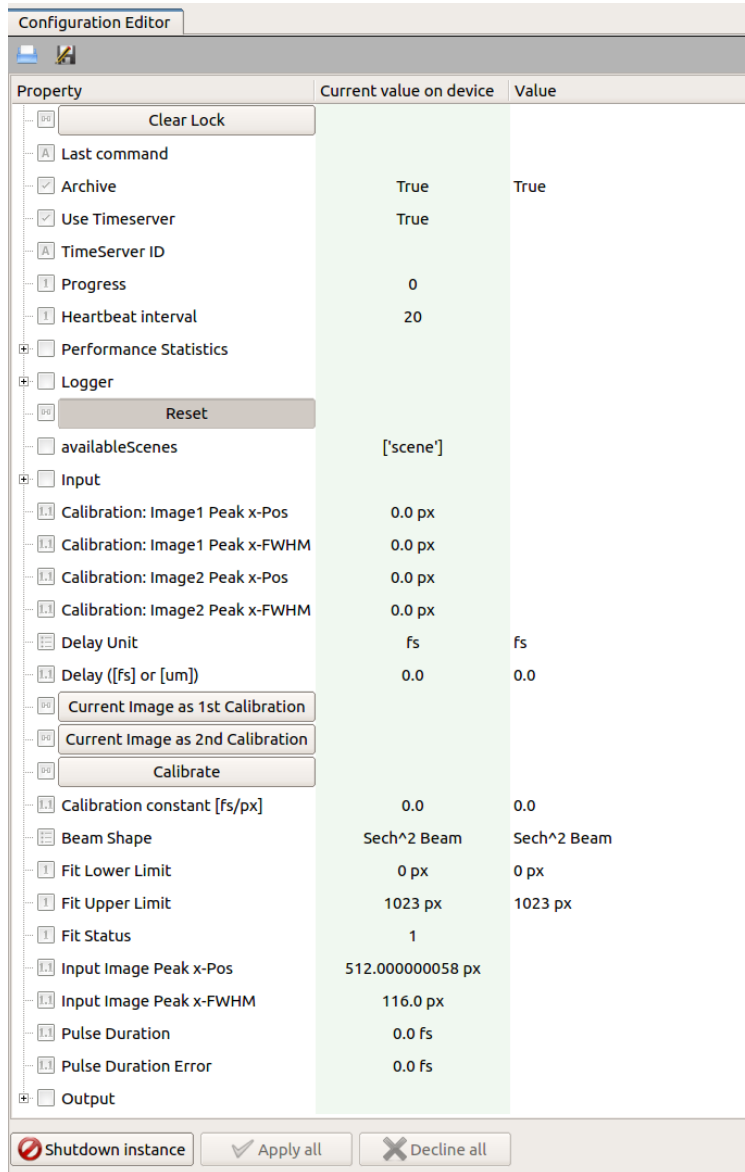


Fig. 1.3: Configuration Editor of the autocorrelator device.

- 4: The iteration is not making good progress, as measured by the improvement from the last five Jacobian evaluations,
- 5: The iteration is not making good progress, as measured by the improvement from the last ten iterations,
- 'unknown': "An error occurred.

The result of pulse duration is presented only in case of a solution is found, and the fit status value is lower than four.

Calibration

To overcome the difficulty in measuring the incident angle ϕ of the primary beams, the following method is applied.

By shifting the mirror stage in the optical delay line, Fig. 2.1, a delay Δt is added between the two input pulses, resulting in a shift ΔZ_0 of the center of SH transverse distribution

$$\Delta Z_0 = \frac{\Delta t \cdot u}{2 \cdot \sin(\phi/2)}$$

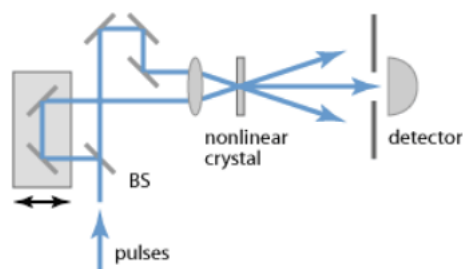


Fig. 2.1: Setup of an intensity autocorrelator. BS refers to the beam splitter.

Combining equations on transverse profile D_z with shift ΔZ_0 the dependence on the intersection angle ϕ is removed, and the pulse duration can be obtained as

$$\tau_p = D_z \cdot \frac{1}{2} \cdot \frac{\Delta t}{\Delta Z_0}$$

The ratio $K = \frac{\Delta t}{\Delta Z}$ is a calibration factor which allows the conversion of the SH transverse profile (measured in pixel units) to the pulse time profile (measured in femtosecond units).

Its determination with sufficient accuracy is challenging. To overcome this difficulty the following procedure is applied. One of the two optical paths can be varied by pulling or pushing one mirror in the line in a controllable way using a micrometer. A change Δl of the micrometer head position results in a pulse delay of $\Delta t = 2\Delta l/c$ and in the shift ΔZ_0 . Thus, shifting the SH distribution, as measured in the CCD camera, in two extreme opposite positions (1 & 2) of the sensitive area allows the measurements of calibration factor with a lower relative uncertainty as shown in the steps below:

$$\Delta t = 2\Delta l/c$$

$$\Delta t_1 - \Delta t_2 = 2(\Delta l_1 - \Delta l_2)/c$$

Considering the above expression of τ_p ,

$$\Delta t_1 - \Delta t_2 = 2 \cdot \tau_p / D_z (\Delta Z_1 - \Delta Z_2)$$

$$(\Delta l_1 - \Delta l_2)/c = \tau_p / D_z (\Delta Z_1 - \Delta Z_2)$$

resulting in

$$\tau_p = D_z \cdot \frac{1}{2} \cdot \left(\frac{2}{c} \cdot \frac{\Delta l_1 - \Delta l_2}{\Delta Z_1 - \Delta Z_2} \right)$$

This way, the calibration factor $K = \left(\frac{2}{c} \cdot \frac{\Delta l_1 - \Delta l_2}{\Delta Z_1 - \Delta Z_2} \right) [ps]$ can be calculated with a larger relative precision using a reproducible and controllable procedure.

It should be noted that the multiplying factor 1/2 in the above equation results from the initial and non-realistic assumption of a rectangular time profile and uniform transverse intensity profile for the incoming beams. More realistic models for the unknown time shape of initial pulses should be considered. Assuming the Gaussian and hyperbolic secant shapes for the pulse time-profile results in the factors 1/2 and 1/1.54, respectively.

The oscillator pulse duration is then calculated as the mean value of these extracted values, and the contribution from model uncertainty to the global systematical uncertainty can be estimated as half of the maximum deviation between the two calculated values.

The above mentioned calibration steps are handled by the device configuration editor. The user should take care to properly select the fitting region reducing the contribution from the fundamental beams. The fitting window can be optimized configuring the keys **Fit Lower Limit** and **Fit Upper Limit**. Also, attention should be taken in order not to cut the profile tail of the SH beam thus affecting the measurement of the FWHM.

After moving the generated SH beam to one side of the sensitive area of the CCD camera (by properly translating the mirror stage in the optical delay line with the micrometer), by clicking on **Current Image as 1st Calibration** the current values of peak position and FWHM will be set as **Image1 Peak (x)** and **Image1 FWHM (x)**, respectively. Similarly, the second set of calibration parameters are obtained steering the SH profile in the other side of the camera and clicking on **Current Image as 2nd Calibration**.

Once the two calibration images are acquired, the calibration constant K can be calculated by clicking on **Calibrate** after setting

- **Delay Unit** to μm ;
- **Delay** to the entire translation of the mirror stage, equivalent to $(\Delta l_1 - \Delta l_2)$. This measurement should be taken by the user;

or, in case the optical delay between the two calibration images was provided already in femtosecond unit, after setting

- **Delay Unit** to fs ;
- **Delay** to the time delay.

The extracted **Calibration constant** allows to calculate the pulse duration from the measured FWHM D_z ,

$$\tau_p = D_z \cdot \alpha \cdot K,$$

α being the multiplication factor originating from the model assumed for the time-profile of the pulse.

The uncertainty of the pulse duration is preliminary estimated via error propagation by the uncertainty on the fit FWHM, assuming the uncertainty of the calibration constant is negligible and that no correlation between the fit parameters exists.

Device Scenes

At the moment, one scene is auto-generated by the device.

It can be opened either by right-clicking on the device name, and selecting from the drop up menu the item *Open device scene*, or double-clicking on the device name.

An example of scene is presented in Fig. 3.1:

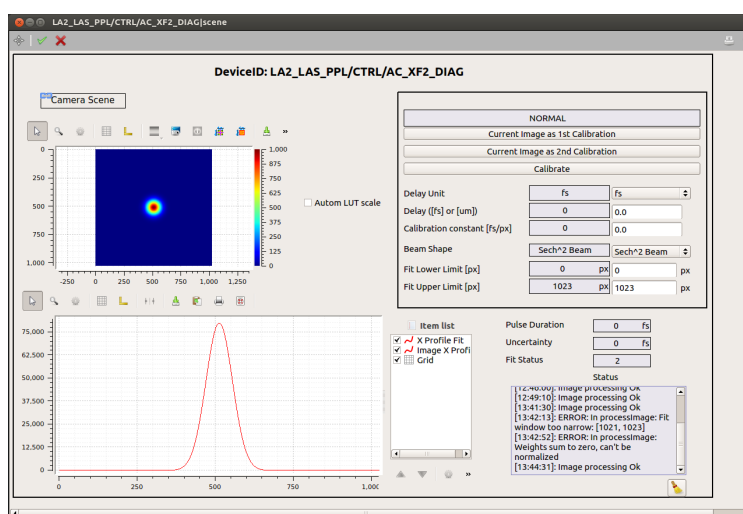


Fig. 3.1: The scene of the auto-correlator device.

All calibration parameters are available in the upper-right sub-panel. The image x-profile is shown superimposed to the fitting function. To deselect one of the graphs use the item list widget. If not yet visible, this widget can be activated from the drop up menu showing up by right-clicking on the graph.

A log of the device **status** is also provided. Note that only messages appeared after the opening of the scene will be displayed.

A link to the camera auto-generated scene is provided, allowing the user to configure the camera without having to navigate in the project.

Troubleshooting

Some typical errors have been identified up to now:

- In case the camera device is not instantiated or it is stopped the peak position and FWHM should be null, and no calculation of the pulse duration can be performed;
- In case no calibration constant is provided, either inserted by the user (if previously known) or by following the calibration procedure described in the text, the pulse duration is not calculated;
- In case the calibration constant is inserted by the user, and the results appear to be very different from what expected, the value used might describe no more the current optical setup of the autocorrelator device. A new calibration measurement could be performed;
- In case the uncertainty arising from the fit procedure is relative large, likely the model used in the fit is not appropriate:
 - try to use a different available model;
 - try to optimize the fitting region;
 - verify that the tails of the second harmonic beam are well within the fitting area;
- In case no available model describes correctly the data, an optimization of the optical line setup could be attempted.

Indices and tables

- `genindex`
- `modindex`
- `search`