

Zyla sCMOS

Version 1.11 rev 10 June 2016



Hardware Guide

	PAGE
SECTION 1: INTRODUCTION.....	8
1.1 HELP AND TECHNICAL SUPPORT	9
1.2 DISCLAIMER.....	10
1.3 TRADEMARKS AND PATENT INFORMATION	10
1.4 SPECIFICATIONS	10
1.5 COMPONENTS	11
1.5.1 ACCESSORIES	11
1.6 POWER AND SIGNAL CONNECTIONS	12
1.6.1 POWER SUPPLY UNIT (PSU)	12
1.6.2 CONNECTORS.....	13
1.6.3 MULTI I/O TIMING CABLE PIN OUTS	13
1.6.4 IMPEDANCE INFORMATION	14
1.6.5 OTHER CONNECTIONS	14
1.7 COOLING HOSE CONNECTORS (WATER COOLED MODELS ONLY)	15
1.7.1 IMPORTANT CONSIDERATIONS WHEN USING COOLING SYSTEMS	15
SECTION 2: FEATURES AND FUNCTIONALITY.....	16
2.1 SCMOS STRUCTURE AND OPERATION.....	16
2.2 ROLLING AND GLOBAL SHUTTER	17
2.2.1 ROLLING SHUTTER.....	18
2.2.2 GLOBAL SHUTTER (Zyla 5.5 ONLY)	19
2.2.3 SELECTING ROLLING OR GLOBAL SHUTTER.....	20
2.2.3.1 EXAMPLES OF TYPICAL APPLICATIONS FOR GLOBAL SHUTTER MODE.....	20
2.2.4 ROLLING AND GLOBAL SHUTTER MECHANISMS	21
2.3 UNDERSTANDING READ NOISE IN sCMOS	22
2.3.1 SPURIOUS NOISE FILTER	23
2.3.2 BLEMISH CORRECTION.....	23
2.4 DUAL AMPLIFIER DYNAMIC RANGE	24
2.5 SENSOR READOUT OPTIMIZATION	26
2.5.1 GAIN CHANNEL CONTROL	26
2.5.2 PIXEL READOUT RATE	27
2.5.3 ROI SUB-IMAGE SETTINGS.....	27

2.6	TRIGGER MODES.....	28
2.6.1	EXAMPLE SYSTEM USAGE OF ACQUISITION EVENTS	29
2.6.2	ROLLING SHUTTER TRIGGERING MODES	30
2.6.2.1	TIMING PARAMETERS AND EXT TRIGGERING FOR ZYLA 5.5	31
2.6.2.2	TIMING PARAMETERS AND EXT TRIGGERING FOR ZYLA 4.2	31
2.6.2.3	ROLLING SHUTTER INTERNAL TRIGGERING (NON-OVERLAP MODE)	32
2.6.2.4	ROLLING SHUTTER INTERNAL TRIGGERING (OVERLAP MODE)	33
2.6.2.5	ROLLING SHUTTER EXTERNAL / SOFTWARE TRIGGERING (NON OVERLAP MODE).....	35
2.6.2.6	ROLLING SHUTTER EXTERNAL EXPOSURE TRIGGERING (NON-OVERLAP MODE)	36
2.6.2.7	ROLLING SHUTTER EXTERNAL EXPOSURE TRIGGERING (OVERLAP MODE).....	37
2.6.2.8	ROLLING SHUTTER EXTERNAL START TRIGGERING	38
2.6.2.9	ROLLING SHUTTER GLOBAL CLEAR INTERNAL (NON-OVERLAP MODE)- ZYLA 4.2 ONLY.....	39
2.6.2.10	ROLLING SHUTTER GLOBAL CLEAR EXTERNAL/SOFTWARE TRIGGERING (NON-OVERLAP MODE) - ZYLA 4.2 ONLY	40
2.6.2.11	ROLLING SHUTTER GLOBAL CLEAR EXTERNAL EXPOSURE TRIGGERING (NON-OVERLAP MODE) - ZYLA 4.2 ONLY	41
2.6.2.12	ROLLING SHUTTER TRIGGERING CONSTRAINTS	42
2.6.3	GLOBAL SHUTTER TRIGGERING MODES (ZYLA 5.5 ONLY).....	43
2.6.3.1	GLOBAL SHUTTER INTERNAL TRIGGERING (NON-OVERLAP MODE)	44
2.6.3.2	GLOBAL SHUTTER INTERNAL TRIGGERING (OVERLAP MODE)	46
2.6.3.3	GLOBAL SHUTTER EXTERNAL/SOFTWARE TRIGGERING	47
2.6.3.4	GLOBAL SHUTTER EXTERNAL EXPOSURE TRIGGERING (NON-OVERLAP MODE).....	49
2.6.3.5	GLOBAL SHUTTER EXTERNAL EXPOSURE TRIGGERING (OVERLAP MODE)	50
2.6.3.6	GLOBAL SHUTTER EXTERNAL START TRIGGERING	51
2.6.3.7	GLOBAL SHUTTER TRIGGERING CONSTRAINTS	52
2.7	ACQUISITION MODES	53
2.7.1	SINGLE SCAN.....	53
2.7.2	KINETIC SERIES.....	53
2.7.3	ACCUMULATE	53
2.7.4	RUN TILL ABORT ACQUISITION.....	53

2.7.4.1	LIVE MODE.....	54
2.7.5	FAST EXPOSURE SWITCH	54
2.7.6	FRAME RATE CONTROL	54
SECTION 3: INSTALLATION		55
3.1	SAFETY CONSIDERATIONS.....	55
3.2	CAMERA MOUNTING.....	55
3.3	COOLANT HOSE CONNECTION AND DISCONNECTION (WATER COOLED MODELS)	56
3.3.1	COOLANT RECOMMENDATIONS.....	56
3.3.2	CONNECTING THE COOLANT HOSES	56
3.3.3	DISCONNECTING THE COOLANT HOSES	56
3.4	RECOMMENDED COMPUTER REQUIREMENTS.....	57
3.5	INSTALLING SOFTWARE (SOLIS OR SDK3)	58
3.6	INSTALLING THE CAMERA FRAMEGRABBER CARD	59
3.7	INSTALLING THE PCIe CARD DEVICE DRIVER (CAMERA LINK MODELS)	60
3.7.1	IMPORTANT INSTALLATION INFORMATION FOR USB 3.0 CAMERAS	60
SECTION 4: OPERATION		61
4.1	CONNECTING YOUR CAMERA AND SWITCHING IT ON.....	61
4.2	USING THE LENS MOUNT ADAPTORS	62
4.2.1	C-MOUNT ADAPTOR.....	62
4.2.2	CS-MOUNT ADAPTOR.....	63
4.2.3	F-MOUNT ADAPTOR	63
4.3	COOLING	64
4.3.1	INTEGRAL THERMAL PROTECTION.....	64
4.3.2	TE COOLING (AIR-COOLED MODELS)	64
4.3.3	TE COOLING (WATER COOLED MODELS)	64
4.4	SOFTWARE SUPPORT.....	65

SECTION 5: MAINTENANCE	66
5.1 REGULAR CHECKS	66
5.2 ANNUAL ELECTRICAL SAFETY CHECKS	66
5.3 FUSE REPLACEMENT	66
5.4 COOLING HOSES AND CONNECTIONS	66
SECTION 6: TROUBLESHOOTING	67
6.1 CAMERA BUZZER DOES NOT SOUND ON START-UP	67
6.2 CAMERA IS NOT RECOGNIZED BY PC	67
6.3 BUZZER SOUNDS CONTINUOUSLY	67
6.4 FAN NOT OPERATING AS EXPECTED	67
6.5 CAMERA DOES NOT COOL TO THE REQUIRED TEMPERATURE	67
6.6 FIRE, AUX_OUT AND ARM OUTPUTS NOT FUNCTIONING CORRECTLY	68
6.7 EXTERNAL TRIGGER INPUT NOT FUNCTIONING CORRECTLY	68
6.8 PREVENTING CONDENSATION	68
APPENDIX A: MECHANICAL DRAWINGS	69
APPENDIX B: DEW POINT GRAPH	71
APPENDIX C: OTHER INFORMATION	72

SAFETY AND WARNING INFORMATION



PLEASE READ THIS INFORMATION FIRST BEFORE USING YOUR ZYLA sCMOS CAMERA.

1. To ensure correct and safe operation of this product, please read this guide before use and keep it in a safe place for future reference
2. If equipment is used in a manner not specified by Andor, the protection provided by the equipment may be impaired
3. Before using the system, please follow and adhere to all warnings, safety, manual handling and operating instructions located either on the product or in this Hardware Manual
4. The Andor Zyla is a precision scientific instrument containing fragile components. Always handle with care
5. Do not expose the product to extreme hot or cold temperatures
6. For air-cooled models, ensure that a minimum clearance of approximately 100mm (4") is maintained in front of all ventilation slots and the fan inlet
7. For water-cooled models, ensure that cooling water supply is connected prior to powering the camera.
8. Do not expose the product to open flames
9. Do not allow objects to fall on the product
10. Do not expose the product to moisture, wet or spill liquids on the product. Do not store or place liquids on the product. If a spillage occurs on the product, switch off power immediately, and wipe off with dry, lint-free cloth. If any ingress has occurred or is suspected, unplug mains cable, do not use, and contact Andor service
11. The product contains components that are extremely sensitive to static electricity and radiated electromagnetic fields, and therefore should not be used, or stored, close to EMI/RFI generators, electrostatic field generators, electromagnetic or radioactive devices, or other similar sources of high energy fields
12. Operation of the system close to intense pulsed sources (e.g. plasma sources, arc welders, radio frequency generators, X-ray instruments, and pulsed discharge optical sources) may compromise performance if shielding of the Zyla is inadequate
13. This product is not designed to provide protection from ionising radiation. Any customer using this product in such an application should provide their own protection
14. The Andor Zyla is for use in research laboratories and other controlled scientific environments
15. This equipment has not been designed and manufactured for the medical diagnosis of patients
16. Use only the power supply cord provided with the system for this unit. Should this not be correct for your geographical area contact your local Andor representative
17. Only the correctly specified mains supply and fuse must be used
18. Make sure the electrical cord is located so that it will not be subject to damage
19. There are no user-serviceable parts beyond the specified user accessible areas of the product and the enclosure must not be opened. Only authorised service personnel may service this equipment



20. The above label indicates that this unit contains components which are sensitive to and can be damaged by electrostatic discharge. When working on a unit which is not enclosed it is necessary to follow anti static precautions to ensure damage does not occur.

REGULATORY COMPLIANCE



The Zyla 4.2 and Zyla 5.5s CMOS Cameras comply with the requirements of the EU EMC and LV Directives through testing to EN 61326-1 and EN 61010-1.



This product requires a DC power supply (refer to **Section 1.6**)

This product complies with international EMC standards, but please note that this includes a permissible loss of performance as allowed by EN 61326 1:2013 and IEC 61326 1:2012 for a defect that rarely occurs during 4 kV contact ESD strikes. This issue manifests itself in the worst case as one row experiencing 5 times the typical noise that persists for the duration of a kinetic series, but disappears thereafter. This issue is theoretical and has never been seen during normal use

REVISION HISTORY

Version	Released	Description
1.0	17 Jul 2012	Initial Release
1.1	29 Oct 2012	General minor updates throughout.
1.2	15 Oct 2013	General enhancements to presentation (all Sections) Text revised to improve clarity of information (all Sections). Zyla Components revised to show standard supplied components more clearly (Section 1.5) Additional accessories revised to match current options available (Section 1.5.1) Additional information added for use for lens mount adaptors (Section 4.2) Updated timing diagrams and tables for Rolling and Global Shutter Modes (Section 2) Updated to show 12-bit data range and gain channels options (Sections 2.6 and 2.7)
1.3	16 Jan 2014	Added Information to support the Zyla 4.2 model (all Sections). Added Rolling Shutter Global Clear modes (Section 2.6) Timing data updated to align with current data (Section 2.6)
1.4	13 Feb 2014	Information added for Water Cooled Zyla models (all Sections). Further information on frame rate performance for each model under different shutter and ROI modes added (Section 2.5)
1.5	09 Jun 2014	Updates to include support for Zyla USB 3.0 model (All Sections) Added new section to cover disabling USB host controller power management (Section 3.7.1) Updated frame rates table for USB 3.0 models (Section 2.5) Updated presentation throughout to match current company branding Updated Figures 4 (Section 1.6.4), 7 (Section 2.1) and 12 (Section 2.4) to enhance quality Included updates to show support for Windows 8 (Sections 3.4 and 3.7)
1.6	15 Sep 2014	Updated accessories list to show that longer lengths of USB 3.0 cable are available, but not supplied or verified by Andor (Section 1.5.1)
1.7	03 Feb 2015	Updated information for Zyla 5.5 model: low vibration fan (as used in Zyla 4.2) fitted to air-cooled units from February 2015 onwards, max. Ambient temperature changed to 30°C. Edited Mechanical drawing for water cooled models to show distance from C-mount to camera faceplate when installed is 6mm nominal.
1.8	17 Nov 2015	Revised Rolling Shutter Sections (Section 2.6.2.3 to 2.6.3.12) Corrected cross-reference to Mechanical Drawings (Section 3.2) Mechanical drawings updated (Appendix A)
1.9	01 Feb 2016	Updated PCI slot for USB Card to x4 or greater- previously was x4 only (Sections 3.4 and 3.6).
1.10	19 May 2016	Added BIOS options checking and setting for Camera Link models used with non-Andor supplied PCs (Section 3.8).
1.11	10 Jun 2016	Added note regarding a potential loss of performance in an ESD strike event (Safety and Warning Section)

SECTION 1: INTRODUCTION

Thank you for choosing the **Zyla Scientific CMOS (sCMOS) camera**. You are now in possession of a revolutionary new **sCMOS camera**, a breakthrough technology based on the next-generation CMOS image sensor (CIS) design and fabrication techniques. The camera offers two image sensors of 5.5 Megapixels (Zyla 5.5) and 4.2 Megapixels (Zyla 4.2). It is available in both air cooled, and water cooled forms and with Camera Link and USB 3.0 connection options.



Figure 1: The Zyla sCMOS Camera. Left: Water Cooled model, right: air cooled model

The Zyla sCMOS camera offers:

- Extremely low noise
- Rapid frame rates
- Wide dynamic range
- High quantum efficiency (QE)
- High resolution
- Large field of view

This manual contains useful information and advice to ensure you get the optimum performance from your new system. If you have any questions regarding your Zyla, camera please feel free to contact Andor directly, or via your local representative or supplier.

1.1 HELP AND TECHNICAL SUPPORT

If you have any questions regarding the use of this equipment, please contact the representative* from whom your system was purchased, or:

Europe

Andor Technology

7 Millennium Way
Springvale Business Park
Belfast
BT12 7AL
Northern Ireland
Tel. +44 (0) 28 9023 7126
Fax. +44 (0) 28 9031 0792

USA

Andor Technology

425 Sullivan Avenue
Suite # 3
South Windsor
CT 06074
USA
Tel. +1 (860) 290 9211
Fax. +1 (860) 290 9566

Asia-Pacific

Andor Technology (Japan)

IS Building
3-32-42 Higashi-Shinagawa
Shinagawa-ku,
Tokyo 140-0002
Japan
Tel: +81 3 6732 8968
Fax: +81 3 6732 8939

China

Andor Technology

Unit 1, Building A,
66 Zhufang Rd, Haidian Dist.
Beijing 100085
China
Tel: +86 (0)10 8271 9066
Fax. +86(0)10 8271 9055

***NOTE:** the contact details for your nearest representative can be found on our website:

<http://www.andor.com/ContactSupport.aspx?type=s>

Introduction

1.2 DISCLAIMER

THE INFORMATION CONTAINED HEREIN IS PROVIDED "AS IS" WITHOUT WARRANTY, CONDITION OR REPRESENTATION OF ANY KIND, EITHER EXPRESS, IMPLIED, STATUTORY OR OTHERWISE, INCLUDING BUT NOT LIMITED TO, ANY WARRANTY OF MERCHANTABILITY, NON-INFRINGEMENT OR FITNESS FOR A PARTICULAR PURPOSE.

IN NO EVENT SHALL ANDOR BE LIABLE FOR ANY LOSS OR DAMAGE, WHETHER DIRECT, INDIRECT, SPECIAL, INCIDENTAL, CONSEQUENTIAL OR OTHERWISE HOWSOEVER CAUSED WHETHER ARISING IN CONTRACT TORT OR OTHERWISE, ARISING OUT OF OR IN CONNECTION WITH THE USE OF THE INFORMATION PROVIDED HEREIN.

COPYRIGHT AND PROTECTIVE NOTICES:

The copyright in this document and the associated drawings are the property of Andor Technology and all rights are reserved. This document and the associated drawings are issued on condition that they are not copied, reprinted or reproduced, nor their contents disclosed.

The publication of information in this documentation does not imply freedom from any patent or proprietary right of Andor Technology or any third party.

1.3 TRADEMARKS AND PATENT INFORMATION

Andor, the Andor logo, Zyla and Solis are trademarks of Andor Technology. Andor is an Oxford Instruments company. All other marks are property of their owners. Changes are periodically made to the product and these will be incorporated into new editions of the manual. New releases of the manual are available through MyAndor: <http://my.andor.com/login.aspx>.

1.4 SPECIFICATIONS

Table 1: Technical Specifications of the Zyla sCMOS Camera

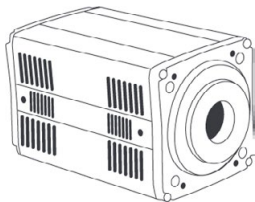
PARAMETER	SPECIFICATION	
AC/DC Power supply ratings	100 - 240 VAC, 43 - 67 Hz	
AC/DC Input Current	1.4 A Max	
AC/DC Output Power	12V DC +/- 5 % @ 5.0A Max (Peak < 6A for < 30 sec after power up)	
Camera Power Consumption	12V @ 5A Max, 12V @ 2.5A Nominal	
Location to be used	Indoor use only	
Altitude	Up to 2000 m	
Ambient Operating Temperature range	Zyla 5.5: 0°C to 30°C (air cooled models*) 0°C to 35°C (water cooled models)	Zyla 4.2: 0°C to 27°C (air-cooled) 0°C to 35°C (water cooled)
Storage temperature	-10°C to +50°C	
Operating relative humidity	< 70% non-condensing	
Overvoltage category	CAT II AC/DC, CAT I Camera	
Pollution degree	Pollution degree 2. Normally only non-conductive pollution occurs. Occasionally, however, a temporary conductivity caused by condensation must be expected.	
Ingress protection rating	IP20	
Electromagnetic compatibility	This is a Class A product. In a domestic environment this product may cause electromagnetic interference, in which case the user may be required to take adequate measures	
Cooling vent clearance	100 mm minimum	
Dimensions	140 x 82 x 80 mm	
Weight	1 kg [2 lb 3 oz]	

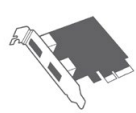





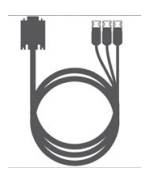


Specifications are subject to change without notice.* Max ambient for Zyla 5.5 aircooled models manufactured prior to Feb 2015 is 35°C.

1.5 COMPONENTS

The standard components supplied with the Zyla are shown in **Table 2**:

Table 2: Standard Components supplied with the Zyla

Description		Quantity
	Zyla 5.5 or Zyla 4.2 sCMOS Camera with Integral Lens Mount Adaptor (C, CS or F-mount: as selected at time of ordering)	1

Description		Quantity	Description		Quantity
	PCIe Frame Grabber Card	1		Anti-static Strap	1
	Camera to PC Interface cables (s) USB 3.0 and 3-tap 10-tap	1 x 3 m 2 x 3 m		Allen Key (1.5 mm)	1
	Power Supply Unit (PS12) and Country specific Power Cord	1		Hardware Guide on CD	1
	Multi I/O Timing cable	1 x 3 m		Quick Start Guide	1
	Coolant hose inserts (for water cooled models only)	2			

1.5.1 ACCESSORIES

There is a range of optional and additional accessories available for your Zyla camera including:

- Mounting Adapters (C-mount, CS-mount or F-mount)
- Re-circulator and Compact Chiller Units (water cooled models must be water cooled via some method)
- Camera Link Cables (5 meter and 10 meter active including power supply)
- Fibre-optic Extenders (30 and 100 meter fibre-optic extender solutions)
- Software Development Kit (SDK)
- Solis Image Capture and Analysis Software
- iQ Imaging Software

Note: Longer lengths of USB 3.0 cable are available from other suppliers such as Stemmer/CEI.

Please contact Andor or your nearest Andor representative for further information.

1.6 POWER AND SIGNAL CONNECTIONS

1.6.1 POWER SUPPLY UNIT (PSU)

The Zyla camera is powered by an external 12 V DC Power Supply Unit (PSU). The Zyla camera is connected via a 4 pin power connector - refer to **Section 1.6.5**. The external power supply has an IEC male socket that requires a certified mains lead with an IEC female plug for connection to the mains electrical supply.

The Zyla PSU is for use with Telecommunications, Computer, Industrial Controller, and OA Systems and must only be used indoors

 The Zyla requires a Direct Current (DC) supply.



Figure 2: Power Supply Unit

NOTES:

1. The electrical mains lead should be certified for use in your country and in applicable countries the plug must be fitted with a 240V 5A fuse.
2. If users use any other power supply, they do so at their own risk.

1.6.2 CONNECTORS



Figure 3: Zyla Back Plate Connections (Left) Cameralink versions (Right) USB 3.0 version

1.6.3 MULTI I/O TIMING CABLE PIN OUTS

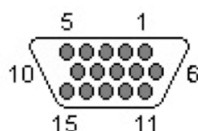


Table 3: Multi I/O Timing Cable Pinouts 15-way D type connector

1	ARM	9	Reserved
2	AUX_OUT_1	10	Reserved
3	FIRE n	11	Reserved
4	FIRE	12	Reserved
5	AUX_OUT_2	13	Reserved
6	Ground	14	Reserved
7	External Trigger	15	Reserved
8	Spare (I)		

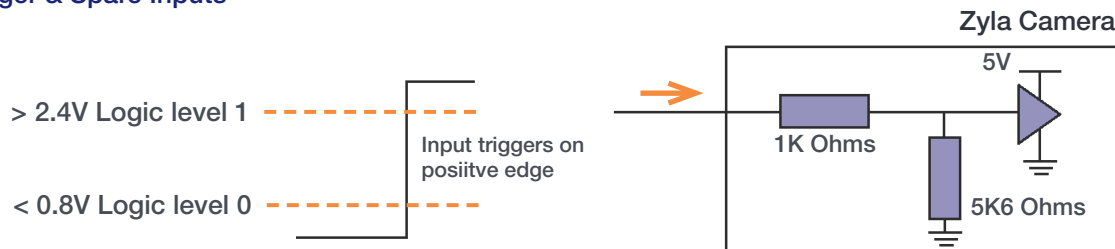
- **External Trigger** and **Spare** input are 5V TTL input. By default they trigger on a rising edge.
- **Fire, Fire n, Arm, AUX_OUT_1 and AUX_OUT_2** outputs are all TTL timing outputs (please also refer to **Section 1.6.4** for information on impedance matching)
- **TTL I/O** can be individually inverted via software (e.g. Solis or SDK)
- **Pins 9 to 15** are reserved and should not be used.
- **AUX_OUT_1** supplies the 'FIRE ALL' output by default. This is the logical AND of the FIRE pulses associated with Row #1 and Row #n (the last row read out in the image frame). Therefore the **FIRE ALL pulse represents the time within a frame when all rows on the sensor are simultaneously exposing**. **AUX_OUT_1** is also configurable as FIRE, FIRE n and FIRE ANY. The **FIRE ANY pulse represents the time within a frame when any row of the image frame is exposing**. Refer to **Section 2.6** for the behaviour of these signals and to the SDK3 manual for configuring the AUX_OUT_1 output.

NOTE: This configurable output is only available on cameras with FPGA version numbers \geq 20120802 and Solis versions \geq 4.22.30001.0 (SDK users require version \geq 3.5.30001.0).

- **AUX_OUT_2** output is reserved for future use.
- I/O Timing Interface cable (Andor part number ACC-ACZ-05612) gives access to all of the above I/O functions (excluding Ground and Reserved).

1.6.4 IMPEDANCE INFORMATION

Ext Trigger & Spare Inputs



Arm, Fire, Fire n and AUX_OUT Outputs

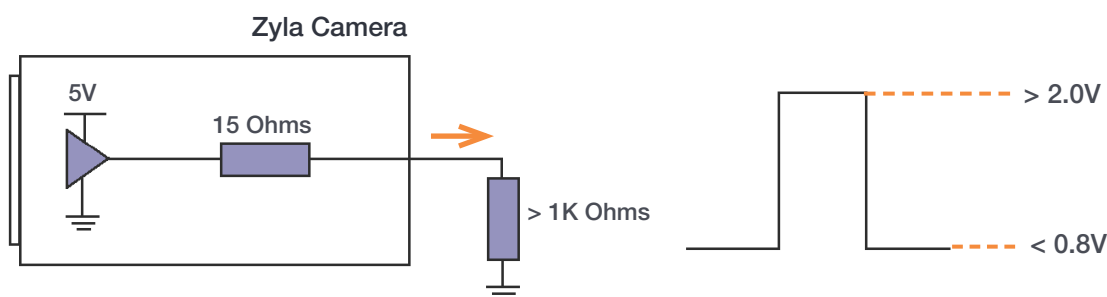


Figure 4: Zyla Connection Impedance Information

1.6.5 OTHER CONNECTIONS

- **CL1:** This is the connection for the 3-tap (single cable) Camera Link version of the Zyla camera.
- **CL2:** This connection is required for the full 10-tap (double cable) Camera Link version of the Zyla 5.5 and Zyla 4.2 camera and is not present on the Zyla 5.5 3-tap version.
- **USB 3.0:** This is the connection for the USB 3.0 version of the Zyla camera. It is not present on the Camera Link versions.
- **Power:** A 4 pin power connector is fitted for power connection, details are shown below:

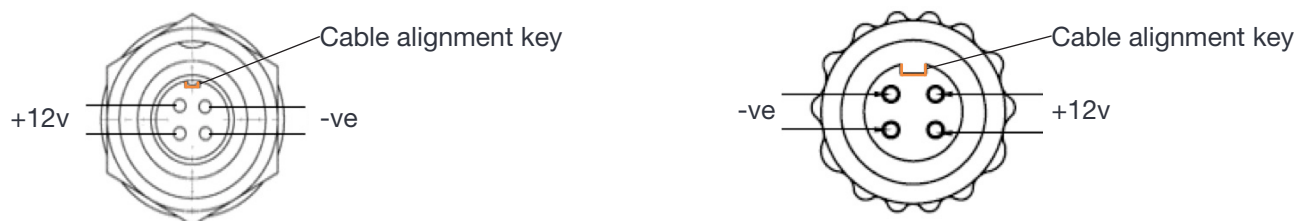


Figure 5: (Left, Camera Connection) and (Right, Connector socket)

WARNING: Ensure that the power switch is in the 'OFF' position and the connector orientation is correct before inserting the power connector. The power connector is 'keyed' - Never forcibly insert the connector otherwise damage to the equipment may occur.

1.7 COOLING HOSE CONNECTORS (WATER COOLED MODELS ONLY)



There are two connectors to allow connection of Water Cooled Zyla models to a water cooler or re-circulator. Hose inserts are provided to enable connection to coolant hoses.



Figure 6: Zyla Cooling Hose Connectors

Please refer to **Section 3.3** for details of connector and hose type compatibility and for connection and disconnection information.

1.7.1 IMPORTANT CONSIDERATIONS WHEN USING COOLING SYSTEMS

- Always ensure that the temperature of the liquid coolant circulated through the camera head is above the dew point of the camera ambient temperature and humidity conditions. Refer to the dew point graph in **Appendix B** for guidance.
- Use of coolant at or below the dew point can result in permanent damage to the camera head, due to formation of condensation on internal components.
- In the event that replacement hose inserts / barbs are required, please contact your local Andor representative.

SECTION 2: FEATURES AND FUNCTIONALITY

2.1 sCMOS STRUCTURE AND OPERATION

sCMOS technology has been developed specifically to overcome many of the limitations that have marred other scientific detector technologies, resulting in an imaging detector that provides exceptional performance for many applications.

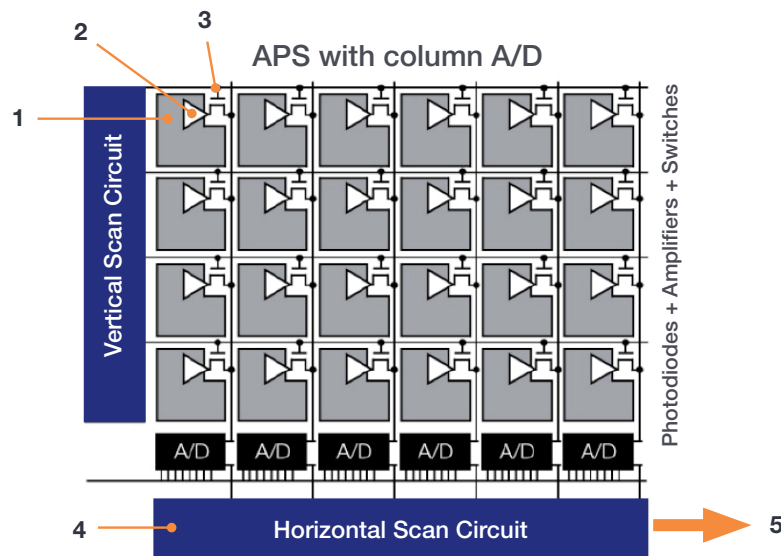


Figure 7: sCMOS Sensor Architecture

As illustrated above, the CMOS sensor is an “Active Pixel Sensor” (APS) whereby each pixel has its own integral amplifier and the sequence of operation is as follows:

1. Light hits sensor and generates charge
2. The photo-generated charge is converted to an analog voltage inside each pixel amplifier
3. Pixel voltage is transferred to the column bus via a row select signal
4. The analog voltage is then converted to a digital signal via columns of A/D (analog to digital) converters.
5. The final digitized signals are then read out sequentially at a pixel readout speed of up to 280 MHz for the Zyla 5.5 and 270 MHz for the Zyla 4.2 (in x2 halves).

NOTES:

The diagram, above is representative- the light sensitive area is contiguous as the photodiodes for each pixel are buried within the sensor. Each Pixel also has a microlens to maximize sensitivity to light.

For Rolling Shutter mode operation, pixels in each row are exposed and the charge converted to a voltage simultaneously before being digitized then read out sequentially

For Global Shutter mode, each pixel in the sensor begins an exposure simultaneously and then ends this exposure simultaneously

2.2 ROLLING AND GLOBAL SHUTTER

The sCMOS sensor used by the Zyla 5.5 offers a choice of both Rolling and Global shutter, providing superior application flexibility. Rolling and Global shutter modes describe two distinct sequences through which the image may be read off a sCMOS sensor. In rolling shutter, charge is transferred from each row in sequence during readout, whereas in global shutter mode each pixel in the sensor effectively ends the exposure simultaneously. However, lowest noise and fastest frame rates are achieved from rolling shutter mode.

Traditionally, most CMOS sensors offer either one or the other, but very rarely does the user have the choice of both from the same sensor. For the Zyla 5.5, it is possible to select between either readout mode from the same sensor, so the most appropriate mode can be selected for the specific application. The Zyla 4.2 offers Rolling Shutter only.

2.2.1 ROLLING SHUTTER

In Rolling Shutter mode, adjacent rows of the array are exposed at slightly different times as the readout ‘waves’ sweep through each half of the sensor. Therefore, each row will start and end its exposure slightly offset in time from its neighbour. In the case of the Zyla 5.5, at the maximum readout rate of 560 MHz (as each half of the sensor is at 280 MHz), this offset between adjacent row exposures is $\sim 10 \mu\text{s}$. The rolling shutter readout mechanism is illustrated in **Figure 8** below. From the point of view of readout, the sensor is split in half horizontally. Rows are read out from the centre outwards, row after row. At the start of an exposure, the wave sweeps through each half of the sensor, switching each row in turn from a ‘keep clean state’, in which all charge is drained from the pixels, to an ‘exposing state’, in which light induced charge is collected in each pixel. At the end of the exposure, the readout wave again sweeps through the sensor, transferring the charge from each row into the readout node of each pixel. The important point is that each row will have been subject to exactly the same exposure time, but the row at the top (or bottom) edge of the sensor would have started and ended its exposure $\sim 10 \text{ ms}$ ($1080 \text{ rows} \times 10 \mu\text{s}/\text{row}$) after the rows at the centre of the sensor (when using 560 MHz readout rate).



Figure 8: Rolling Shutter Exposure and Readout

Rolling shutter can be operated in a ‘continuous’ mode when capturing a kinetic series of images, whereby after each row has been read out it immediately enters its next exposure. This ensures a 100% duty cycle, meaning that no time is wasted between exposures and, perhaps more importantly, no photons are wasted. At the maximum frame rate for a given readout speed (e.g. 100 fps at 560 MHz for the Zyla 5.5) the sensor is continuously reading out, i.e. as soon as the readout fronts reach the top and bottom of the sensor, they immediately return to the centre to readout the next exposure.

A potential downside of rolling shutter is spatial distortion resulting from the above described exposure mechanism. This has historically been more apparent in devices such as CMOS camcorders, where the entire image field could be moved (for example by the user rapidly panning the camera) at a rate that the image readout could not match; thus, objects could appear at an angle compared to their actual orientation. In reality, despite the time-offset readout pattern, rolling shutter mode is appropriate for the majority of scientific applications, especially where the exposure time is equal to or greater than the sensor readout time, discussed later.

2.2.2 GLOBAL SHUTTER (ZYLA 5.5 ONLY)

Global shutter mode, which can also be thought of as a ‘snapshot’ exposure mode, means that all pixels of the array are exposed simultaneously. In most respects, global shutter can be thought of as behaving like an Interline CCD sensor. Before the exposure begins, all pixels in the array will be held in a ‘keep clean state’, during which charge is drained into the anti-bloom structure of each pixel. At the start of the exposure each pixel simultaneously begins to collect charge and is allowed to do so for the duration of the exposure time. At the end of exposure each pixel transfers charge simultaneously to its readout node. Importantly, global shutter can be configured to operate in a continuous ‘overlap’ mode (analogous to Interline CCD), whereby an exposure can proceed while the previous exposure is being readout out from the readout nodes of each pixel. In this mode, the sensor has a 100% duty cycle, again resulting in optimal time resolution and photon collection efficiency.

However, the mechanism of global shutter mode demands that a reference readout is performed ‘behind the scenes’, in addition to the actual readout of charge from each pixel. Due to this additional reference readout, global shutter mode carries the trade-off of halving the maximum frame rate that would otherwise have been achieved in rolling shutter mode. In addition, global shutter also increases the RMS read noise by a factor of 1.41 over rolling shutter readout.

Figure 9 below shows a simplified illustration showing sequence of events in global shutter mode:

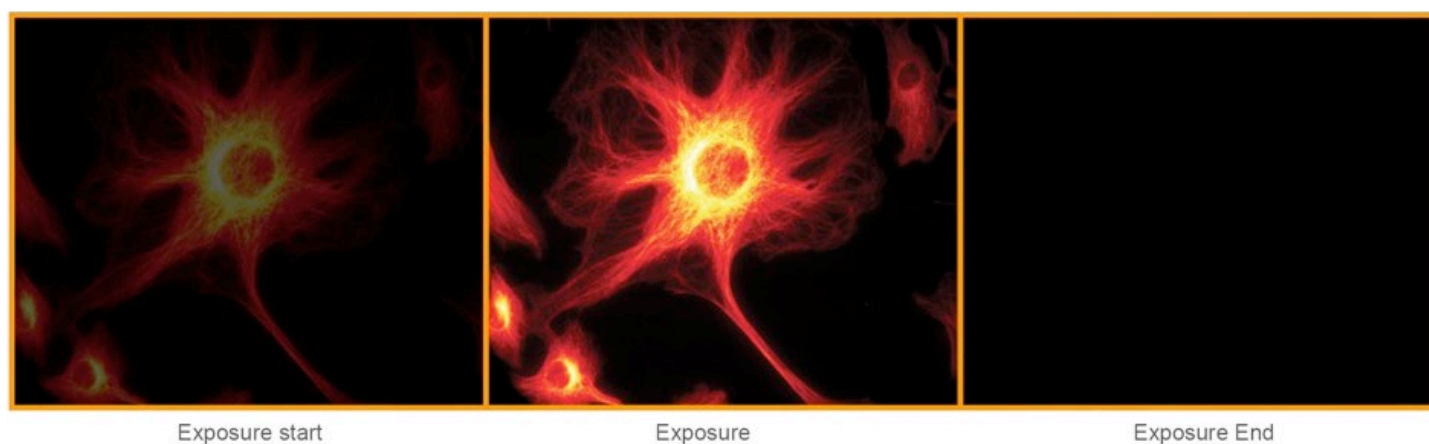


Figure 9: Global Shutter Exposure and Readout

2.2.3 SELECTING ROLLING OR GLOBAL SHUTTER

The selection of Rolling Shutter or Global Shutter modes for the Zyla 5.5 depends on your specific experimental conditions. A summary of the key parameters for each mode is shown in **Table 4**.

Table 4: A Comparison of Rolling and Global Shutter Modes

PARAMETER	ROLLING SHUTTER MODE	GLOBAL SHUTTER MODE (ZYLA 5.5 ONLY)
Frame Rate	Maximum available	Maximum frame rate is halved
Read Noise	Lowest	Increased by 1.41
Spatial Distortion	Dependent on object dynamics and frame rate	None

Rolling Shutter Mode: with the enhanced frame rates and lower noise, is likely to suit the majority of scientific applications. As long as the frame rate is such that the camera is temporally oversampling object dynamics within the image area, negligible spatial distortion will be observed. Such oversampling is good imaging practice, since it is undesirable to have an object travel a significant distance during a single exposure.

Global Shutter Mode: for some specific applications global shutter will be viewed as a necessity. These are shown in **Section 2.2.3.1**.

Refer also to Andor Technical Note, “Rolling and Global Shutter”.

2.2.3.1 EXAMPLES OF TYPICAL APPLICATIONS FOR GLOBAL SHUTTER MODE

- **Applications that require ‘microsecond’ time gating synced to a pulsed light source:** e.g. Laser Induced Breakdown Spectroscopy (LIBS). Global readout involves a step that simultaneously transfers the signal charge of each pixel into the corresponding readout node for that pixel. This transfer step is 2 μ s (at 560 MHz readout rate), facilitating fast exposure end, i.e. ‘Electronic gating’
- **‘Double Exposure’ applications:** e.g. Particle Imaging Velocimetry (PIV), which requires that two back-to-back exposures are acquired with minimal time separation between them. The global shutter 2 μ s transfer time into the readout node defines the minimum time between two consecutive exposures
- **Applications that require exact time correlation between two (or more) points of an image that are separated vertically within the image:** In rolling shutter it takes 10 μ s per row for the 2x readout fronts to move across the image from the centre outwards, reading out one row at a time. At 560 MHz pixel readout rate, this represents 10 ms to cover the distance from centre to outermost rows. That means an object at the centre of the image will begin and end an exposure ~10 ms before an object located at the very top or bottom (although remember that each object will be subject to the same overall exposure time). If a particular application requires that ‘moving or changing’ objects separated by relatively large distances (vertically) be subject to the same beginning and end of exposure, then global shutter mode is required
- **Applications where the entire field of view is fast moving (relative to exposure time):** e.g. high speed machine vision inspection, such as PCB inspection

2.2.4 ROLLING AND GLOBAL SHUTTER MECHANISMS

In Rolling Shutter mode, charge transfer happens on a per row basis whilst in global shutter charge transfer happens for the whole sensor or globally. To read out a pixel in Rolling Shutter mode, the following occurs within the analog circuitry:

1. The read out node is reset
2. The node level (reference level) is measured
3. Charge is transferred from photodiode to node
4. The node level (signal level) is measured
5. Reference level (step 2) is subtracted from signal level (Step 4) to get real signal

This process is commonly referred to as CDS (Correlated Double Sampling) and is done in the analog domain before digitization. The reason it is required is due to what is known as reset noise, this arises because every time the node is reset it does not settle at exactly the same level and hence the actual level must be measured (**Step 2.**) and subtracted from the signal level (**Step 4**) to get the real signal.

Rolling Shutter Mode: charge transfer happens on a per row basis; therefore each row follows steps 1 – 5, until the entire sensor is read out. The disadvantage of this is that the start and end exposure time moves by the row read out time for each subsequent row. So whilst each row of pixels is exposed for exactly the same length of time they do not all start and end at exactly the same time.

Global Shutter Mode: the start and end of the exposure do occur at exactly the same time for every pixel (not just for pixels in the same row); therefore **Step 3.** has to occur for all the pixels at the same time. Because of this, the reset and reference read occur before this global transfer for every row. Since the same read out circuitry is used for every row there is nowhere to store the measured reference level for every pixel and so a reference frame is actually digitized and read out from the sensor and then the signal is digitized and read out from the sensor. The two are subtracted to get the 'real signal'. Reading two frames to get a real signal frame effectively halves the Cycle Time when compared to Rolling Shutter.

2.3 UNDERSTANDING READ NOISE IN sCMOS

sCMOS technology boasts an ultra-low read noise floor that significantly exceeds that of even the best CCDs, and at several orders of magnitude faster pixel readout speeds. For those more accustomed to dealing with CCDs, it is useful to gain an understanding of the nature of read noise distribution in CMOS imaging sensors.

CCD architecture is such that the charge from each pixel is transferred through a common readout structure, at least in single output port CCDs, where charge is converted to voltage and amplified prior to digitization in the Analog to Digital Converter (ADC) of the camera. This results in each pixel being subject to the same readout noise. However, CMOS technology differs in that each individual pixel possesses its own readout structure for converting charge to voltage. In the sCMOS sensor, each column possesses dual amplifiers and ADCs at both top and bottom (facilitating the split sensor readout). During readout, voltage information from each pixel is fed directly to the appropriate amplifier/ADC, a row of pixels at a time (see Technical Note on Rolling and Global Shutter modes).

As a consequence of each pixel having its own individual readout structure, the overall readout noise in CMOS sensors is described as a distribution, as exemplified in **Figure 10** below, which is a representative noise histogram from a Zyla 5.5 camera at the fastest readout speed of 560 MHz (or 280 MHz x 2 halves). It is standard to describe noise in CMOS technology by citing the median value of the distribution. In the data presented, the median value is 1.38 electron RMS. This means that 50% of pixels have a noise less than 1.38 electrons, and 50% have noise greater than 1.38 electrons. While there will be a small percentage of pixels with noise greater than 2 or 3 electrons, observable as the low level tail towards the higher noise side of the histogram, it must be remembered that a CCD Interline camera reading out at 20 MHz would have 100% of its pixels reading out with read noise typically ranging between 6 and 10 electrons RMS (depending on camera design).

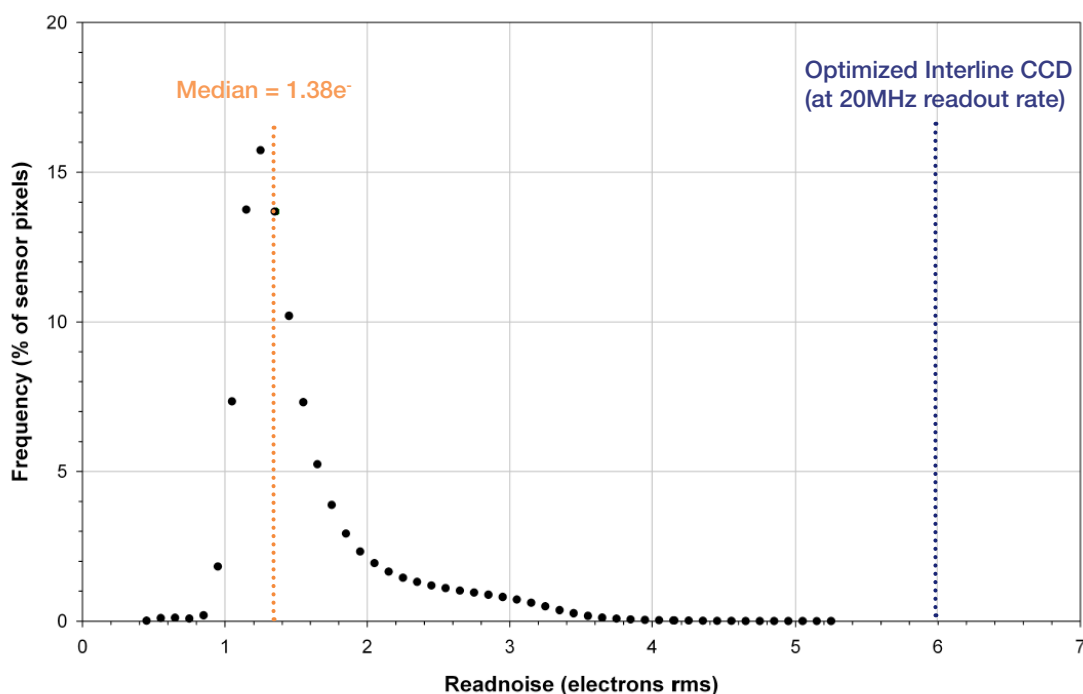


Figure 10: Representative histogram showing read noise distribution at fastest readout speed of Zyla 5.5, 560 MHz (280 MHz x2). The median value of 1.38 e⁻ means 50% pixels have read noise less than 1.38 e⁻ and 50% have greater than 1.38 e⁻. The line at 6 e⁻ represents a typical read noise value from a well optimized Interline CCD – all pixels in a CCD essentially share the same noise value

2.3.1 SPURIOUS NOISE FILTER

The **Spurious Noise** filter corrects for pixels that would otherwise appear as spurious ‘salt and pepper’ noise spikes in the image. The appearance of such noisy pixels is analogous to the situation of Clock Induced Charge (CIC) noise spikes in EMCCD cameras, in that the overall noise of the sensor has been reduced to such a low level, that the remaining small percentage of spurious, high noise pixels can become an aesthetic issue. The filter actively corrects such high noise pixels, replacing them with the mean value of the neighbouring pixels. The filter can be switched on and off by the user prior to data acquisition.

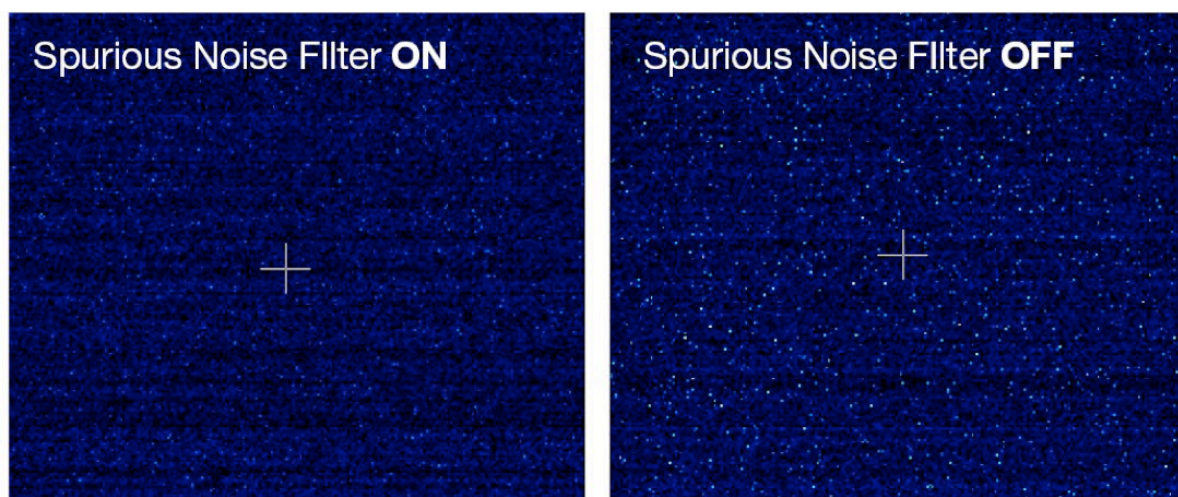


Figure 11: Demonstration of Spurious Noise Filter (Filter On left, Filter Off, Right) on a dark image, 20 ms exposure time, 200 MHz (x2 halves) readout speed

2.3.2 BLEMISH CORRECTION

This **Blemish Correction** filter identifies and compensates for three types of blemishes during the FPGA processing step:

1. Hot Pixel's
2. Noisy Pixel's
3. Unresponsive Pixel's

sCMOS sensors are particularly susceptible to hot pixel blemishes. These are spurious noise pixels that have significantly higher darkcurrent than the average. Through deep TE cooling of the sensor (e.g. -30°C in the Neo), it is possible to dramatically minimize the occurrence of such hot pixels within the sensor, meaning that these pixels can still be used for useful quantitative imaging. However, if deep cooling cannot be achieved it is necessary to use interpolative filters to minimize the hot pixel blemishes. These filters work by taking the mean of the surrounding 8 pixel values and replacing this hot pixel blemish with this mean value. Such interpolation over pixel blemishes can be detrimental in some applications that depend on total quantitative integrity over a limited set of pixels, for example in localization based super-resolution microscopy (such as PALM and STORM techniques) and astronomy. In these applications it is essential for the user to be able to switch off interpolative corrections.

Furthermore, having access to the location of these blemishes allows an accurate map of ‘good’ pixels to be determined by the user. A new service allows the end user to request a ‘hot pixel map’ of their sCMOS sensor from Andor. This map will be generated based on the experimental conditions outlined by the end user.

From the latest general release of Andor SDK3 (version 3.7.30004) and Solis (version 4.24.30004) blemish correction can be switched on and off by the user. Refer to the SDK and Solis User Guide and help information for instructions.

2.4 DUAL AMPLIFIER DYNAMIC RANGE

The Dual Amplifier architecture of the sCMOS sensor in Zyla eliminates the need to choose between low noise or high capacity, in that signal can be sampled simultaneously by both high gain and low gain amplifiers. As such, the lowest noise of the sensor can be harnessed alongside the maximum well depth, affording the widest possible dynamic range. Traditionally, scientific sensors including CCD, EMCCD, ICCD and CMOS, demand that the user must select ‘upfront’ between high or low amplifier gain (i.e. sensitivity) settings, depending on whether they want to optimise for low noise or maximum well depth. Since the true dynamic range of a sensor is determined by the ratio of well depth divided by the noise floor detection limit, then choosing either high or low gain settings will restrict dynamic range by limiting the effective well depth or noise floor, respectively.

For example, consider a large pixel CCD, with 16-bit Analog to Digital Converter (ADC), offering a full well depth of 150,000 e⁻ and lowest read noise floor of 3 e⁻. The gain sensitivity required to give lowest noise is 1 e⁻/ADU (or ‘count’) and the gain sensitivity required to harness the full well depth is 2.3 e⁻/ADU, but with a higher read noise of 5 e⁻. Therefore, it does not automatically follow that the available dynamic range of this sensor is given by $150,000/3 = 50,000:1$. This is because the high sensitivity gain of 1e⁻/ADU that is used to reach 3 e⁻ noise means that the 16-bit ADC will top out at 65,536 e⁻, well short of the 150,000 e⁻ available from the pixel. Therefore, the actual dynamic range available in ‘low noise mode’ is $65,536/3 = 21,843:1$. Conversely, the lower sensitivity gain setting means that the ADC will top out at ~ 150,000 e⁻, but the higher read noise of 5 e⁻ will still limit the dynamic range to $150,000/5 = 30,000:1$ in this ‘high well depth mode’. The sCMOS sensor offers a unique dual amplifier architecture, meaning that signal from each pixel can be sampled simultaneously by both high and low gain amplifiers. The sensor also features a split readout scheme in which the top and bottom halves of the sensor are read out independently. Each column within each half of the sensor is equipped with dual column level amplifiers and dual analog-to-digital converters, represented by the block diagram below:

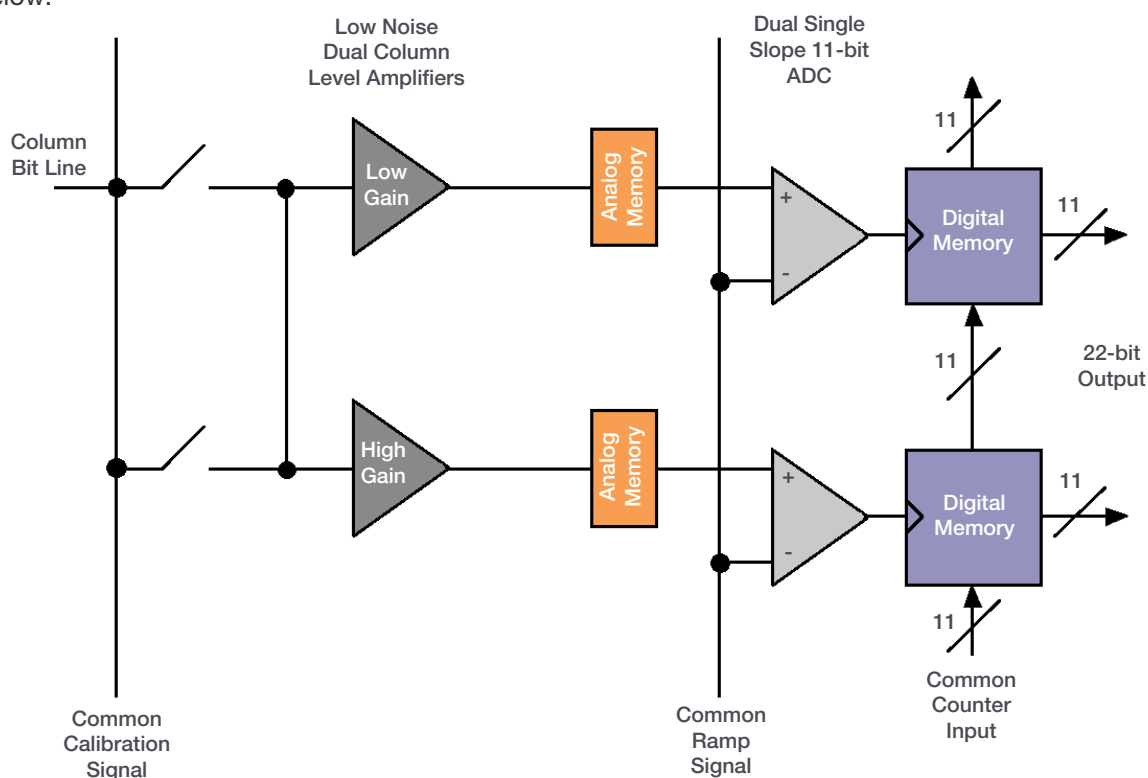


Figure 12: Amplifiers and ADC of the sCMOS Sensor

The dual column level amplifier/ADC pairs have independent gain settings, and the final image (see Figure 13) is reconstructed by combining pixel readings from both the high gain and low gain readout channels to achieve a wide intra-scene dynamic range, uniquely so considering the relatively small 6.5 μm pixel pitch.

The method of combining signals from two 11-bit ADCs can be divided into four basic steps.

1. At the end of the analog chain the “Signal” voltage is applied to two independent amplifiers: the high gain amplifier and the low gain amplifier. This results in two separate digital data streams from the sensor
2. The camera selects which data stream to use on a pixel per pixel, frame by frame basis using a threshold method
3. The data is then compensated for DC offset and gain. Again, this is done on a pixel by pixel basis using the compensation data associated with the data stream. The gain corrects for pixel to pixel relative sensitivity, pixel node amplifier and the high and low amplifier relative gains
4. The pixels are then combined into a single 16-bit image for transfer to the PC

The user maintains the choice of opting to stay with 12-bit single gain channel data if dynamic range is not critical, resulting in smaller file sizes. This in turn offers faster frame rates when continuously spooling through the Camera Link interface and writing to hard disk.

Table 5: Typical performance of supported gain settings of the Zyla 5.5 (Jan 2012 onwards)

AMPLIFIER GAIN (CURRENT ANDOR SDK / SOLIS DESCRIPTION)	MODE	SENSITIVITY e^- / ADU (TYPICAL)	DATA RANGE	EFFECTIVE PIXEL SATURATION LIMIT / e^-	SPOOLING FILE SIZE (PER FRAME)
12-bit (high well capacity)	GS/RS	7.5	12-bit	30,000	8 Mb
12-bit (low noise)	GS	0.42	12-bit	1,700	8 Mb
12-bit (low noise)	RS	0.28	12-bit	1,100	8 Mb
16-bit (low noise and high well capacity)	GS	0.45	16-bit	30,000	10.5 Mb
16-bit (low noise and high well capacity)	RS	0.45	16-bit	30,000	10.5 Mb

Table 6: Typical performance of supported gain settings of the Zyla 4.2

AMPLIFIER GAIN (CURRENT ANDOR SDK / SOLIS DESCRIPTION)	MODE	SENSITIVITY e^- / ADU (TYPICAL)	DATA RANGE	EFFECTIVE PIXEL SATURATION LIMIT / e^-	SPOOLING FILE SIZE (PER FRAME)
12-bit (high well capacity)	RS	7.5	12-bit	30,000	6.1 Mb
12-bit (low noise)	RS	0.28	12-bit	1,100	6.1 Mb
16-bit (low noise and high well capacity)	RS	0.45	16-bit	30,000	8.1 Mb

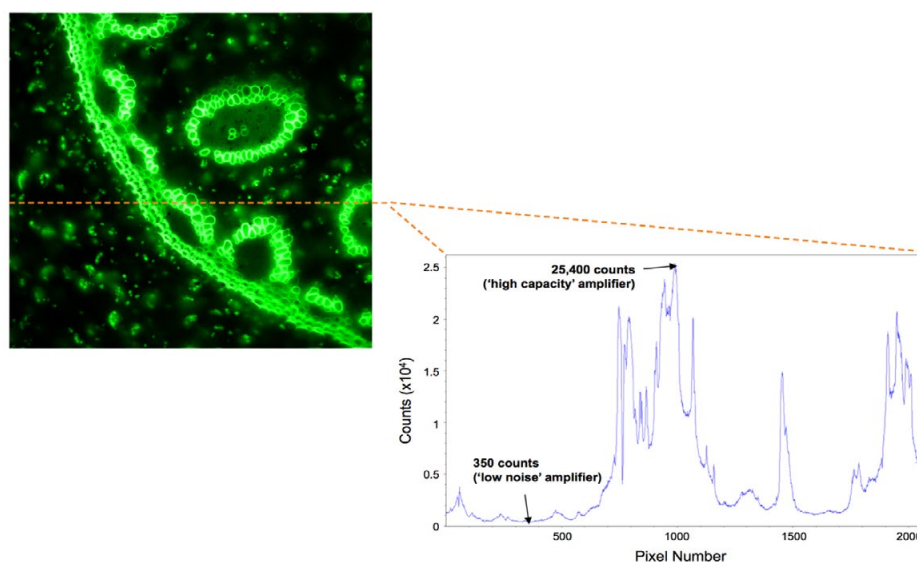


Figure 13: High contrast image of fixed labelled cell. Intensity line profile through single row demonstrates pixel regions that were sampled by high gain (low noise) and low gain (high capacity) amplifiers.

2.5 SENSOR READOUT OPTIMIZATION

To allow the camera to be optimized for the widest range of applications it is important to have flexibility in the readout options available, some of these include the following:

- Gain Channel Control
- Pixel Readout Rate
- ROI sub image settings
- Triggering / Synchronization options (see **Section 2.6**)

The effect of different shutter modes, readout speed and different ROIs on frame rates is shown in **Table 7**.

Table 7: Maximum sustainable frame rates for all electronic shutter modes, readout speeds and gain channel combinations for various (vertically centred) ROIs

Region of Interest	Shutter Mode	Bit-depth	Zyla 5.5 3-tap		Zyla 5.5 10-tap		Zyla 4.2 10-tap		Zyla 5.5 USB 3.0		Zyla 4.2 USB 3.0	
			200 MHz	560 MHz	200 MHz	560 MHz	216 Mhz	540 MHz	200 MHz	560 MHz	216 Mhz	540 MHz
2560 x 2160	Rolling	12-bit	30	30	36	100	-	-	36	40	-	-
		16-bit	22	22	36	75	-	-	30	30	-	-
	Global	12-bit	18	30	18	49	-	-	18	40	-	-
		16-bit	18	22	18	49	-	-	18	30	-	-
2048 x 2048	Rolling	12-bit	38	39	38	105	40	100	38	53	40	53
		16-bit	29	29	38	98	40	100	38	40	40	40
	Global	12-bit	19	39	19	52	-	-	19	52	-	-
		16-bit	19	29	19	52	-	-	19	52	-	-
1920 x 1080	Rolling	12-bit	72	79	72	199	76	192	72	107	76	107
		16-bit	59	59	72	199	76	192	72	80	76	80
	Global	12-bit	35	79	35	97	-	-	35	98	-	-
		16-bit	35	59	35	97	-	-	35	80	-	-
512 x 512	Rolling	12-bit	152	419	152	419	161	403	152	419	161	403
		16-bit	152	419	152	419	161	403	152	419	161	403
	Global	12-bit	73	201	73	201	-	-	73	201	-	-
		16-bit	73	201	73	201	-	-	73	201	-	-
128 x 128	Rolling	12-bit	596	1639	596	1639	631	1578	596	1639	631	1578
		16-bit	596	1639	596	1639	631	1578	596	1639	631	1578
	Global	12-bit	260	716	260	716	-	-	260	716	-	-
		16-bit	260	716	260	716	-	-	260	716	-	-

NOTE: The write speed of the PC hard drive can impose a further restriction on the maximum sustainable frame rate.

2.5.1 GAIN CHANNEL CONTROL

The Zyla offers the user a choice of two 12-bit gain channels (i.e. high or low gain) or a combined '16-bit' setting. The user can choose to stay with 12-bit single gain channel data if dynamic range is not critical, resulting in smaller file sizes. This in turn offers faster frame rates when continuously spooling through the Camera Link interface and writing to hard disk. The 16-bit dual gain channel should be chosen if dynamic range is more important than having the fastest frame rates. Please refer to **Section 2.4** for more detailed explanation of the dual amplifier dynamic range operation.

2.5.2 PIXEL READOUT RATE

The Pixel Readout Rate defines the rate at which pixels are read from the sensor. The faster the readout rate the higher the frame rate that can be achieved. The ability to change the pixel readout rate is important to achieve the maximum flexibility of camera operation.

Slower readout typically allows lower read noise but at the expense of slower frame rates. The following readout rates are available on the Zyla (**Table 8** below shows the typical read noise at each readout rate):

- **Zyla 5.5:** 200 MHz (100 MHz x 2 halves) and 560 MHz (280 MHz x 2 halves)
- **Zyla 4.2:** 216 MHz (108 MHz x 2 halves) and 540 MHz (270 MHz x 2 halves)

Table 8: Typical Read Noise for Rolling and Global Shutter modes

MODEL	RATE	READ NOISE (e ⁻) ROLLING SHUTTER (TYPICAL)	READ NOISE (e ⁻) GLOBAL SHUTTER (TYPICAL)
Zyla 5.5	200 MHz	1.2	2.6
	560 MHz	1.45	2.6
Zyla 4.2	216 MHz	0.9	N/A
	540 MHz	1.1	N/A

Please refer to **Section 2.3** for more information on read noise and the camera performance sheet for read noise values at the various readout speeds.

2.5.3 ROI SUB-IMAGE SETTINGS

ROI Sub Image allows for readout of a particular sub-area of the sensor. When a sub image has been defined, only data from the selected rows will be digitized.

Selecting a sub image increases the frame readout rate and reduces image storage requirements. Examples of sub-image selection and spooling rates are shown in **Table 7**.

It should be noted that for small ROIs (e.g. < 512 x 512 pixels), the achievable frame rate for 3-tap and 10-tap becomes identical. This is due to the fact that the amount of data transmitted across the Camera Link interface for these ROIs is relatively small and the limiting factor becomes the rate at which data can be read off the sensor.

It should also be noted that while ROIs can also be set up for any area of the sensor, those which are vertically centred on the sensor mid line will result in the maximum frame rates.

2.6 TRIGGER MODES

The Zyla camera has the following triggering modes:

- **Internal Trigger** - the camera determines the exact time when an exposure happens based on the acquisition settings entered by the user. This is the most basic trigger mode and requires no external intervention.
- **External Trigger** - the camera and software are in a high state of readiness to accept a trigger from an external source. Refer to **Table 10** for the minimum pulse width required to guarantee a trigger. The external trigger is fed via the External Trigger input on the I/O Connector on the camera head.
- **Software Trigger** works in the a similar manner to External Trigger mode whereby the camera and software are in a high state of readiness and can react extremely quickly to a trigger event issued via software. This mode is particularly useful when the user needs to control other equipment between each exposure and does not know in advance how long such control will take or if the time taken changes randomly.
- **External Start** is a mode where the camera will wait for one external trigger event to occur after the acquisition sequence has been started. Once this external trigger event is detected, the camera will start the Internal Trigger read out process and will progress as if the camera was in internal trigger mode.
- **External Exposure Trigger** is a mode of operation where the exposure time and cycle time are controlled by the external trigger input.

The TTL inputs and outputs may be used to synchronize the camera operation with external events or equipment.

The individual outputs are described in **Sections 2.6.2** and **2.6.3**.

The AUX_OUT_1 output can be configured via software (Solis or SDK) to provide one of the following outputs: FIRE, FIRE n, FIRE ALL, or FIRE ANY.

The default state provides 'FIRE ALL on this output.

The polarity of the TTL inputs and outputs can also be inverted (individually) via either Solis or SDK.

NOTE: 'Row 1' is the first row read out in the image frame. 'Row n' is the last row read out in the image frame.

NOTE: The trigger diagrams in the following sections are for outlining the events and timing of outputs in the various trigger modes and not to scale.

2.6.1 EXAMPLE SYSTEM USAGE OF ACQUISITION EVENTS

Software Acquisition Events are only accessible via SDK- these are not available in Solis, iQ or other software but may be used internally. Refer to the SDK3 manual for further information on configuration of Software Acquisition Events.

Compare the two diagrams below: The first diagram (**Figure 14**) shows the situation prior to the introduction of the software events. In this case the user had to wait until the image frame was completely transferred to the PC before they received any notification that the exposure had completed. This results in a cycle time of approximately 79 ms or a frame rate of 12.6 fps, for this example. With acquisition software events (shown in **Figure 15**) and in particular the ExposureEndEvent the user will be notified as soon as the exposure is complete in advance of readout completion. This in conjunction with faster SW trigger, means that the next acquisition can be started much sooner, resulting in an improvement in frame rate, 25 fps. In addition, the Fire ALL signal can be used to control light sources/z-stages.

Example: 3-tap, RS, 560MHz
30ms exposure, full image

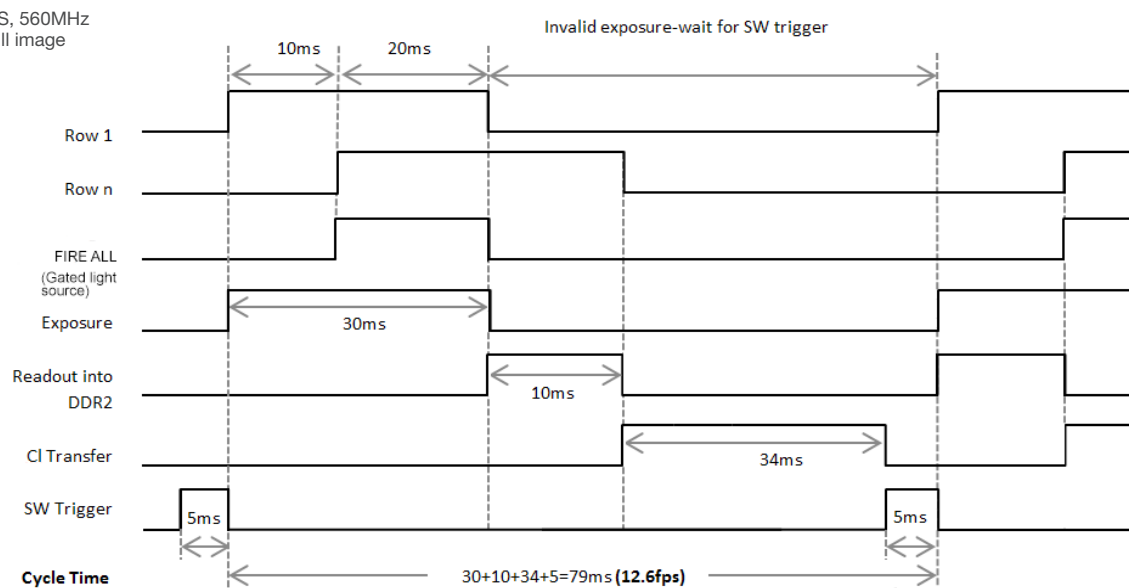


Figure 14: Rolling Shutter - Original Acquisition using SW Trigger

Example: 3-tap, RS, 560MHz
30ms exposure, full image

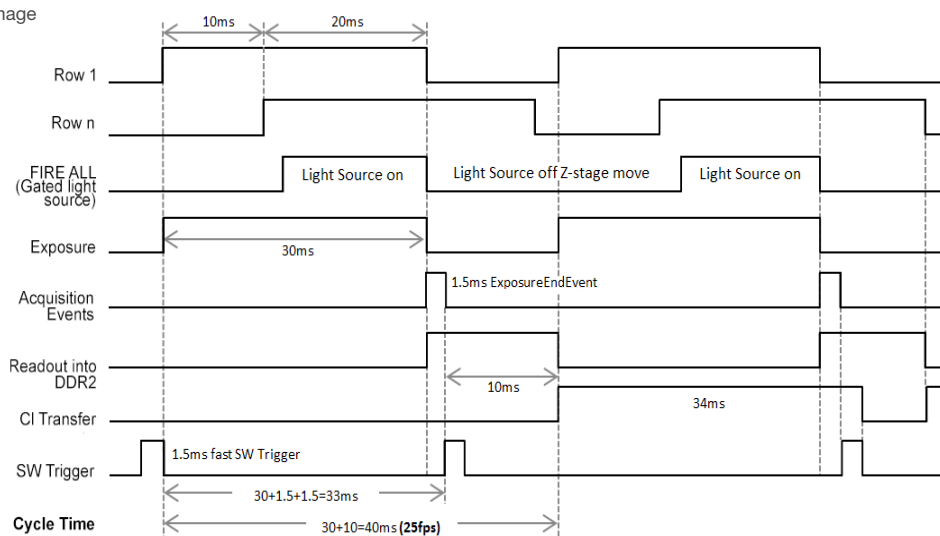


Figure 15: Rolling Shutter - New Acquisition using Events and Faster SW Trigger

2.6.2 ROLLING SHUTTER TRIGGERING MODES

In Rolling Shutter, charge transfer happens on a per row basis:

1. The node is reset,
2. A 'reference' measurement taken, and then the charge transferred,
3. The 'signal' is measured,
4. The reference and signal measurements are then subtracted in the analog domain.
5. Each following row is processed until the whole sensor is read out.

The disadvantage of this is that the start and end exposure time moves by the row read out time for each subsequent row. So whilst each row of pixels is exposed for exactly the same length of time they do not all start and end at exactly the same time.

The Rolling Shutter signals in the diagrams are as follows:

- **Acquisition Start:** This is an internal pulse purely for illustrative purposes and indicates when the camera receives a command from software to start the pre-programmed acquisition sequence
- **FIRE:** (Exposure for Row 1): In Rolling Shutter mode, the FIRE output from the camera indicates to the user the exposure time for the first row
- **FIRE Row n:** (Exposure for Row n): The exposure for Row 2 is delayed by one row time relative to Row 1, Row 3 is delayed by one row time relative to Row 2, etc. for all rows in the frame (up to Row n) - the rows within the selected ROI which are furthest from the sensor centre. This signal is connected to an external output from the camera known as FIRE Row n.
- **FIRE ALL:** The Fire ALL output from the camera indicates when all rows within a frame are being simultaneously exposed.
- **FIRE ANY:** The FIRE ANY output indicates when any row within a frame is being exposed.
- **ARM:** The ARM output from the camera is used in external and software triggering modes to indicate when the camera is ready to accept an incoming trigger. If ARM is low when a trigger event occurs, it will be ignored
- **Global Clear:** The entire sensor can be held in a global clear state which ensures that there is no charge build-up on the sensor
- **Prescan Readout Cycle:** Once an acquisition sequence has been started, the camera is placed into a special cleaning cycle called 'Prescan Readout Cycle' which ensures that charge build up on the sCMOS sensor is kept to a minimum while waiting for the trigger event. The Prescan Readout Cycle consists of repeatedly reading a virtual row from the sensor. Once the trigger is received, a new exposure phase is initiated.
- **Charge Transfer:** This signal indicates when charge in the pixel is transferred to the measurement node and effectively ends the exposure. The charge is transferred while the pulse is HIGH.
NOTE: For clarity this signal is not shown in Rolling Shutter diagrams.
- **Frame Readout Phase:** This signal shows the period during which the signal frame is read out from the sensor
- **EXT Start Delay:** This is the delay between the start of the External Trigger pulse and the start of exposure of row 1.
- * : Marks the start of an exposure.
- * : Marks the end of exposure.

2.6.2.1 TIMING PARAMETERS AND EXT TRIGGERING FOR ZYLA 5.5

The timing tables accompanying each of the triggering diagrams that follow indicate the exposure and cycle times achievable in each triggering mode for the Zyla 5.5. These are based on Frame and Row Periods as shown below.

Table 9: Timing Parameters based on Sensor Clock Speed (Zyla 5.5)

PARAMETER	SENSOR READ OUT RATE	
	200 MHz	560 MHz
1 Row (2624 clock cycles)	25.41 μ s	9.24 μ s
1 Full Frame (2160 rows)	27.44 ms	9.98 ms

The time taken to read out a full row is equal to 2624 clock cycles. The sensor is split into 2 halves with each having an independent data output from the sensor. This means the Frame Period is 1080 rows x 2624 clock cycles. For sub-images the frame readout time is related to the number of rows being displayed.

In External and External Start Triggering Modes, the minimum trigger pulse width detected by the Zyla 5.5 camera is shown below in **Table 10**:

Table 10: Minimum EXT Trigger Width (Zyla 5.5)

PARAMETER	SENSOR READ OUT RATE	
	200 MHz	560 MHz
EXT Trig Pulse Width (2 clock cycles)	19.37 ns	7.04 ns

2.6.2.2 TIMING PARAMETERS AND EXT TRIGGERING FOR ZYLA 4.2

The timing tables accompanying each of the triggering diagrams that follow indicate the exposure and cycle times achievable in each triggering mode for the Zyla 4.2. These are based on Frame and Row Periods as shown below.

Table 11: Timing Parameters based on Sensor Clock Speed (Zyla 4.2)

PARAMETER	SENSOR READ OUT RATE	
	216 MHz	540 MHz
1 Row (2592 clock cycles)	24.00 μ s	9.60 μ s
1 Full Frame (2048 rows)	24.54 ms	9.83 ms

The time taken to read out a full row is equal to 2592 clock cycles. The sensor is split into 2 halves with each having an independent data output from the sensor. This means the Frame Period is 1024 rows x 2592 clock cycles. For sub-images the frame readout time is related to the number of rows being displayed.

In External and External Start Triggering Modes, the minimum trigger pulse width detected by the Zyla 4.2 camera is shown below in **Table 12**:

Table 12: Minimum EXT Trigger Width (Zyla 4.2)

PARAMETER	SENSOR READ OUT RATE	
	216 MHz	540 MHz
EXT Trig Pulse Width (2 clock cycles)	20 ns	7.4 ns

2.6.2.3 ROLLING SHUTTER INTERNAL TRIGGERING (NON-OVERLAP MODE)

Internal Trigger Mode allows the user to configure an exposure time and cycle time. For Internal Triggering Non-overlap mode, the exact acquisition sequence depends on the exposure time and cycle time set as shown in **Figure 16**. The following diagrams show the behaviour of TTL outputs 'Fire', 'Fire n', 'Fire ALL' and 'Fire ANY'.

Fire ALL indicates the time period within a frame during which all rows are exposing simultaneously.

Fire ANY indicates the time period within a frame during which any row is exposing

Initially, the entire sensor is held in a global clear state to ensure that there is no charge build-up on the sensor. Global Clear goes LOW and a frame read out is initiated. This frame is discarded as it does not contain the correct exposure period. Reading out a frame effectively begins a new exposure. When the exposure period has completed, a signal frame read out phase begins. When the frame has been read out completely, the Global Clear is held HIGH until the user-defined cycle time is achieved.

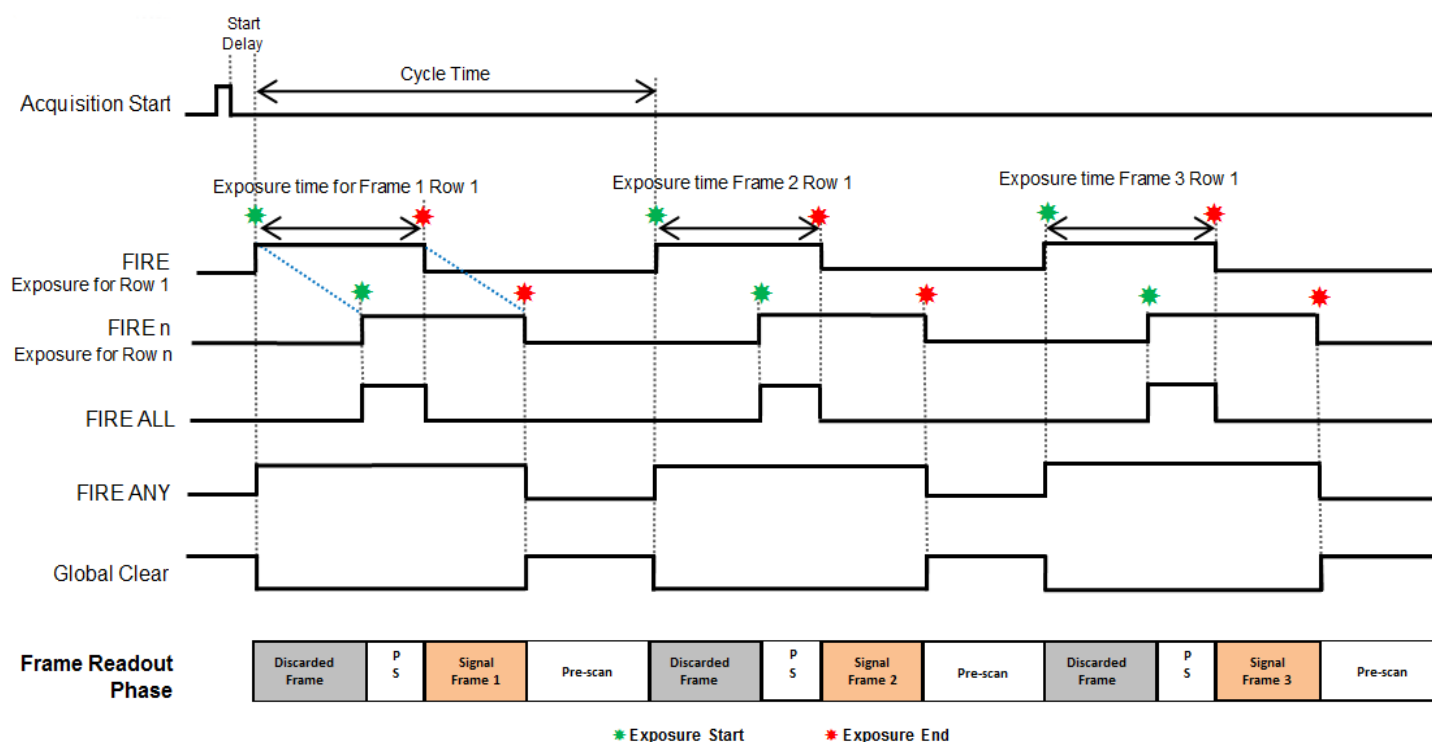


Figure 16: Rolling Shutter Internal Triggering

Table 13: Rolling Shutter Internal Triggering (Long Exposures) Timing Parameters

PARAMETER	MINIMUM	MAXIMUM
Exposure	1 Row	30 s
Cycle Time (1/Frame Rate)	Exposure + 1 Frame + 1 Row	20,000 s
Acquisition Start Delay	-	1 Row

2.6.2.4 ROLLING SHUTTER INTERNAL TRIGGERING (OVERLAP MODE)

Internal Triggering in Overlap Mode allows the user to perform an exposure and acquire images from the sensor simultaneously. This is achieved by starting a new exposure for a new frame while the current frame's exposure is being read out from the sensor. The cycle time is the same as the exposure time entered by the user in this scenario.

Short Exposures

When the required exposure time is less than the time it takes to read out a frame (Short Exposures), the cycle time is always defined by the time taken to read out a frame.

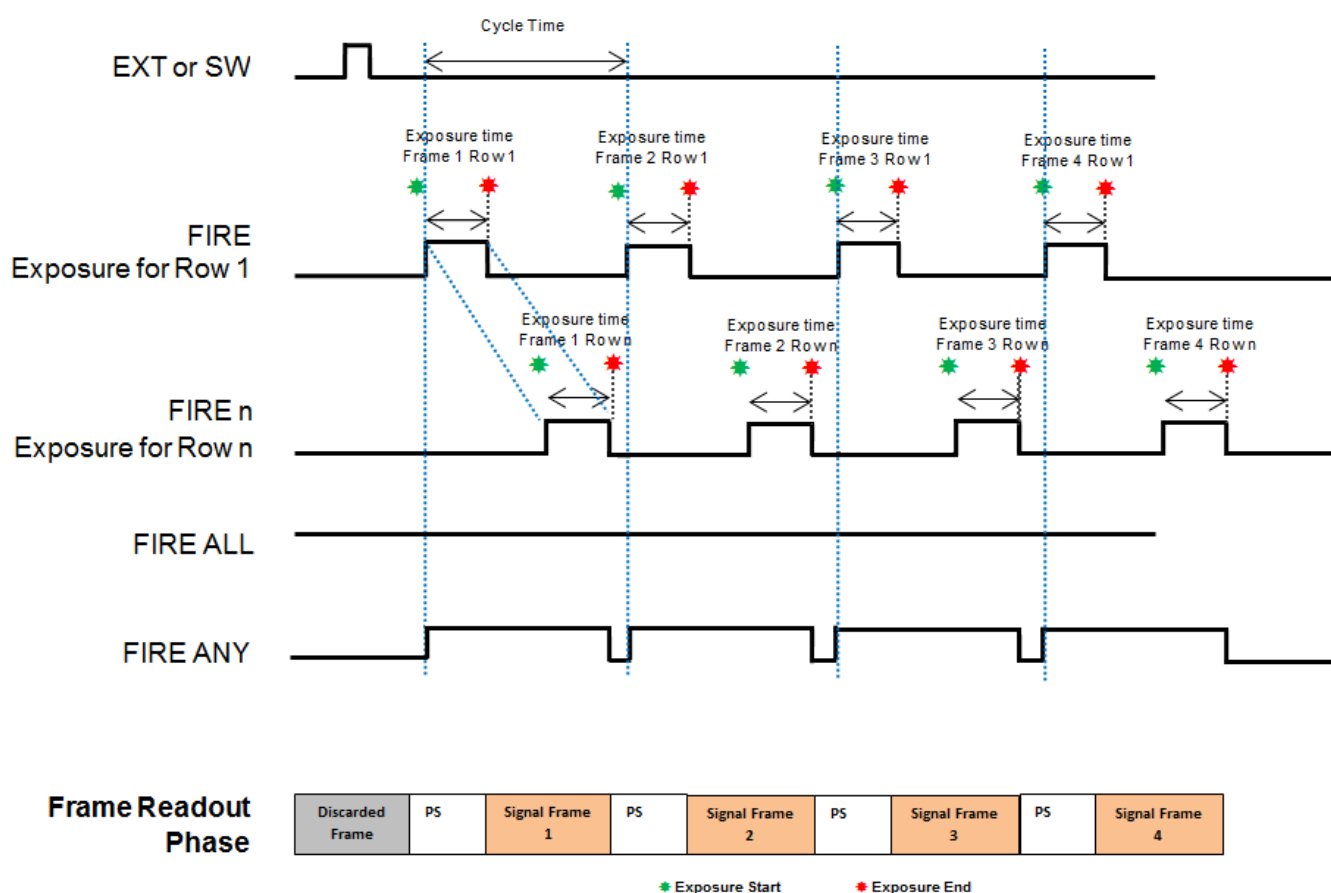


Figure 17: Rolling Shutter Internal Triggering (Short Exposures)

Table 14: Rolling Shutter Internal Triggering (Overlap Mode- Short Exposures) Timing Parameters

PARAMETER	MINIMUM	MAXIMUM
Exposure	1 Row	1 Frame - 1 Row
Cycle Time (1/Frame Rate)	1 Frame	1 Frame + Exposure + 1 Row
FIRE any low period	1 Row	-

Long Exposures

When the required exposure time is greater than the time it takes to read out a frame (Long Exposures), the cycle time is defined by the exposure time.

Initially, the entire sensor is held in a global clear state to ensure that there is no charge build-up on the sensor. Global Clear goes LOW and the first frame read out is initiated. This frame is discarded as it does not contain the correct exposure period. Reading out this first frame effectively begins the first exposure. When the exposure period has completed, a signal frame read out phase begins. As each row is read out, the new exposure for that row begins.

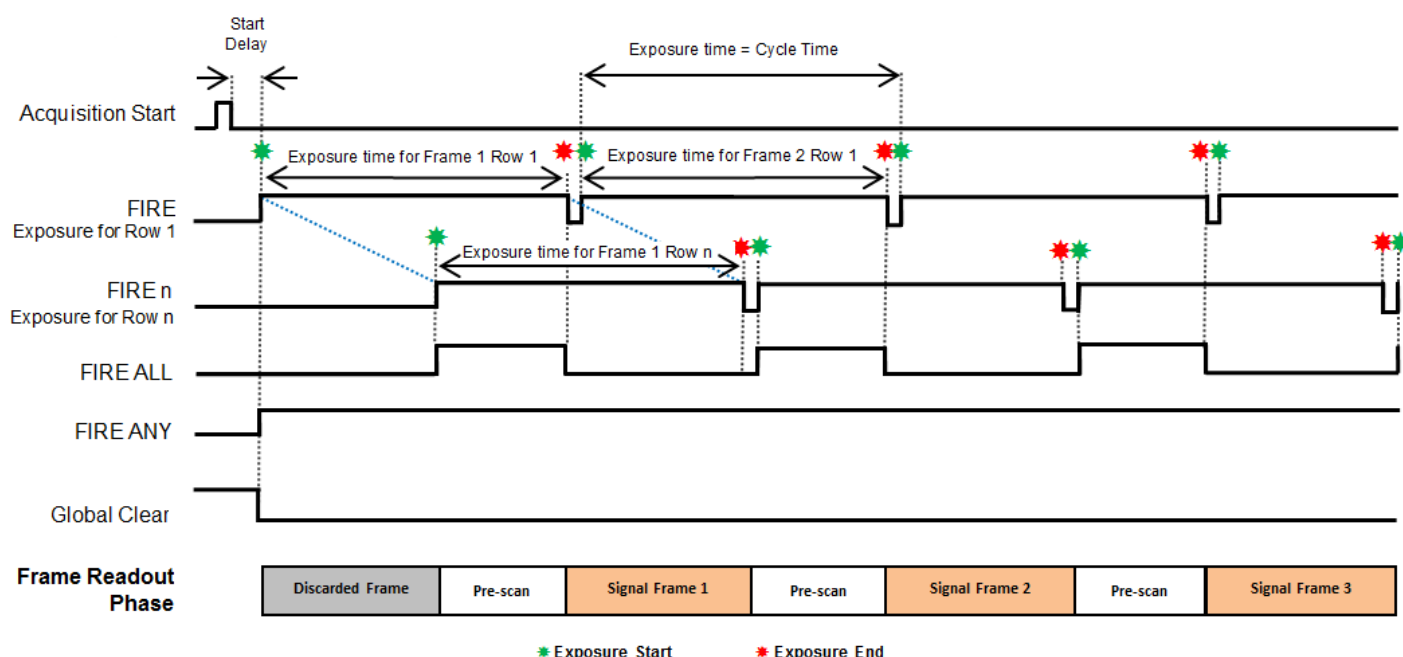


Figure 18: Rolling Shutter Internal Triggering (Overlap Mode) (Long Exposures)

Table 15: Rolling Shutter Internal Triggering (Overlap Mode- Long Exposures) Timing Parameters

PARAMETER	MINIMUM	MAXIMUM
Exposure	1 Row	30 s
Cycle Time (1/Frame Rate)	Exposure Time	Exposure Time + 1 Frame
FIRE low period	1 Row	-

2.6.2.5 ROLLING SHUTTER EXTERNAL / SOFTWARE TRIGGERING (NON OVERLAP MODE)

In this section, both External and Software Trigger are described in the same diagram as the acquisition sequence is the same. The trigger event can either be from the EXT Trigger input or sent via software. While waiting on the trigger event, the sensor is put into a “pre-scan read out cycle” and is held in a Global Clear State which ensures that charge build up on the sensor is kept to a minimum while waiting for the trigger event. The ARM signal is asserted to indicate it is ready to detect an incoming trigger input.

Once the trigger event is detected, Global Clear goes LOW and a frame read out is initiated. This frame is discarded as it does not contain the correct exposure period. Reading out a frame effectively begins a new exposure. When the exposure period has completed, a signal frame read out phase begins. When the frame has been read out completely, the Global Clear is held HIGH until the next trigger event is detected.

The external trigger is fed via the EXT Trigger input on the I/O Connector on the camera head.

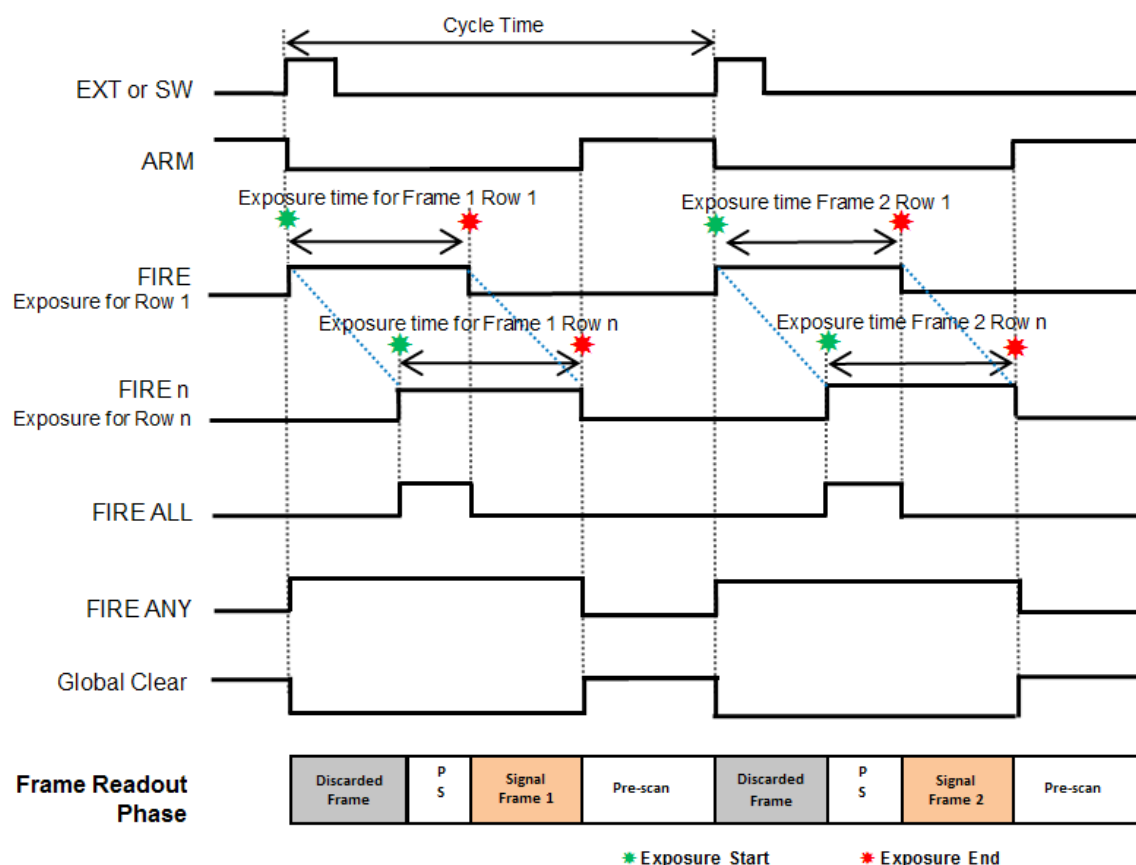


Figure 19: Rolling Shutter External/Software Trigger (Non Overlap Mode)

Table 16: Rolling Shutter External/Software Triggering Timing Parameters

PARAMETER	MINIMUM	MAXIMUM
Exposure	3 Rows	30 s
Cycle Time (1/Frame Rate)	Exposure + 1 Frame + 1 Row	-
External Start Delay	0	1 Row
EXT Trig Pulse Width	2 Sensor Speed Clock Cycles	-

2.6.2.6 ROLLING SHUTTER EXTERNAL EXPOSURE TRIGGERING (NON-OVERLAP MODE)

While waiting on the trigger event, the sensor held in a global clear state and is put into a “pre-scan read out cycle”. On detection of the trigger event, Global Clear goes LOW and a frame read out is initiated. This frame is discarded as it does not contain the correct exposure period. Reading out a frame effectively begins a new exposure. When the external trigger input goes LOW, a signal frame read out phase begins. When the frame has been read out completely, the Global Clear is held HIGH until the next trigger event is detected.

The external trigger pulse width defines the exposure time for all rows, but is only coincident with the exposure time for Row 1. The exposure for Row 2 will be delayed by one row time relative to Row 1 and so forth.

The period of the external trigger pulse defines the overall cycle time. If the width of the trigger event is less than 3 Rows, the falling edge will be missed and a subsequent falling edge will be required to end the exposure.

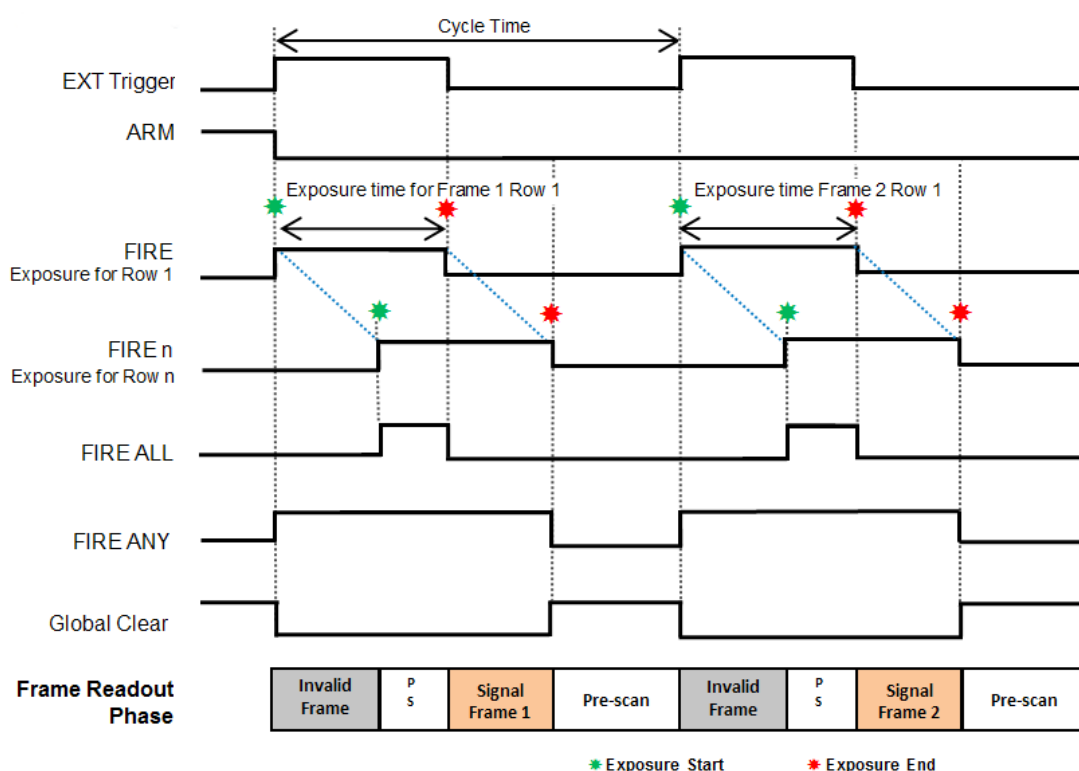


Figure 20: Rolling Shutter External Exposure Triggering (Non Overlap Mode)

Table 17: Rolling Shutter External Exposure Triggering (Non Overlap Mode) Timing Parameters

PARAMETER	MINIMUM	MAXIMUM
Exposure	3 Rows	30 s
Cycle Time (EXT Trig Period)	Exposure + 1 Frame + 1 Row	-
External Start Delay	0	1 Row
EXT Trig Pulse Width	3 Rows	30 s

2.6.2.7 ROLLING SHUTTER EXTERNAL EXPOSURE TRIGGERING (OVERLAP MODE)

In overlap mode, every positive edge of an external trigger will trigger a frame read out and start a new exposure for the next frame. The period of external trigger pulse defines exposure and cycle time for each frame read out.

While waiting on the positive edge of the external trigger, the sensor is held in a global clear state and put into a “pre-scan read out cycle”. On detection of the positive edge, Global Clear goes LOW and a frame read out is initiated. This frame is discarded as it does not contain the correct exposure period. Reading out this first frame effectively begins the first exposure. When the next positive edge of the external trigger is detected, a signal frame read out phase begins. As each row is read out, the new exposure for that row begins.

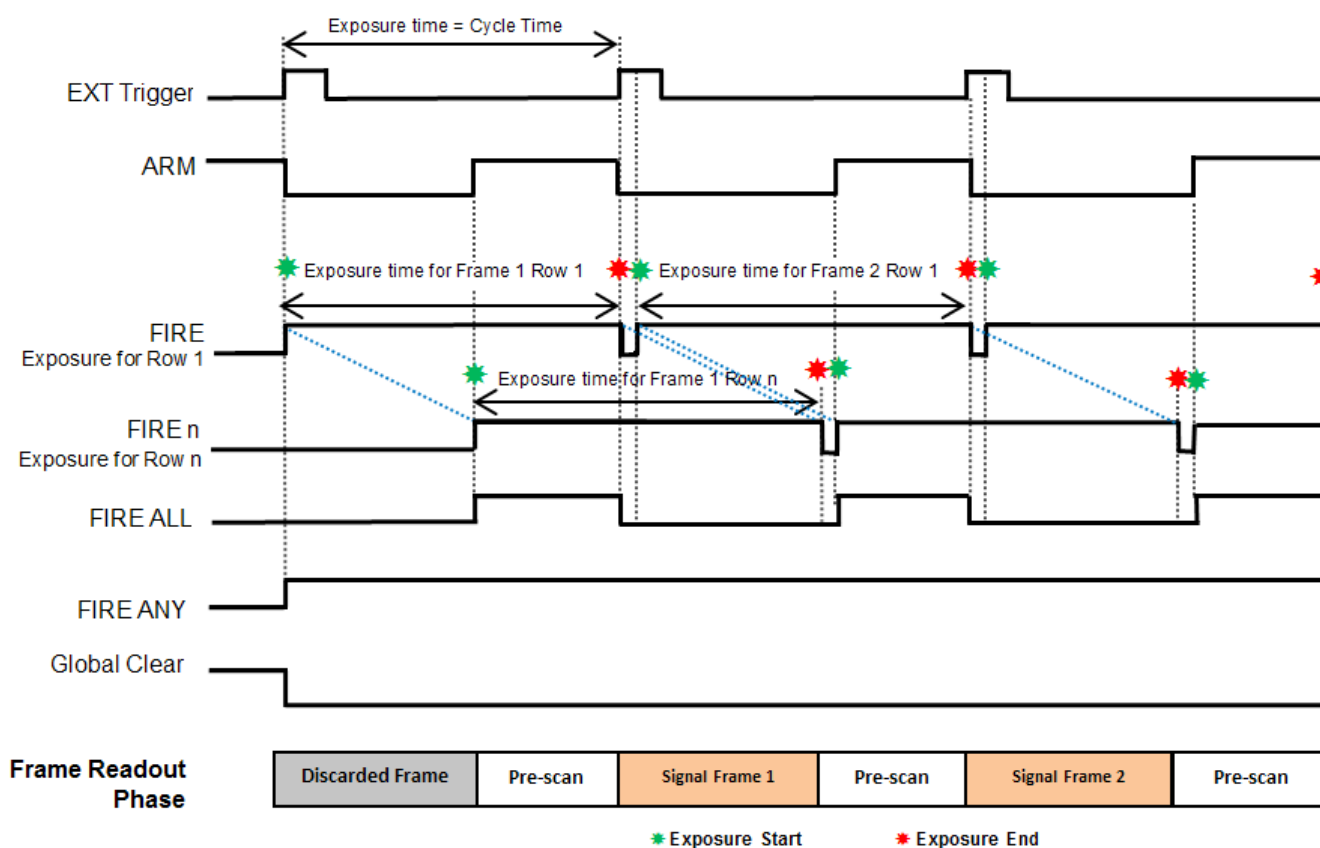


Figure 21: Rolling Shutter External Exposure Triggering (Overlap Mode) (Long Exposures)

Table 18: Rolling Shutter External Exposure Triggering (Overlap Mode) Timing Parameters

PARAMETER	MINIMUM	MAXIMUM
Exposure	1 Frame + 1 Row	30 s
Cycle Time (EXT Trig Period)	Exposure	Exposure
External Start Delay	0	1 Row
EXT Trig Pulse Width	2 Sensor Speed Clock Cycles	-
FIRE low period	20 Sensor Speed Clock Cycles	-

2.6.2.8 ROLLING SHUTTER EXTERNAL START TRIGGERING

In this mode the camera will wait for a single external trigger event. Once this external trigger event is detected, the camera will progress as if the camera was in internal trigger mode (see Section 2.6.2.3 and 2.6.2.4). The ARM signal indicates to the user when the camera is ready to detect an External Start Trigger. Figure 23 below shows the External Start used in Non-overlap Mode, Long Exposure. The delay from the External Trigger to start of exposure is 1 Row.

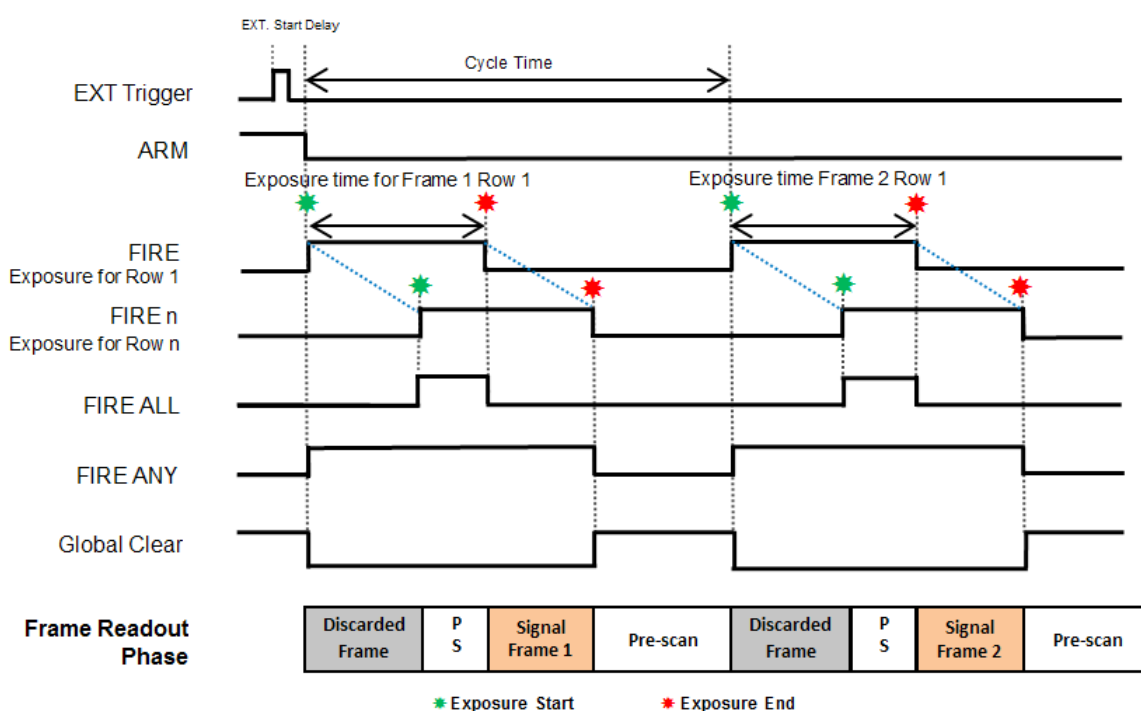


Figure 22: Rolling Shutter External Start Triggering (Non Overlap)

2.6.2.9 ROLLING SHUTTER GLOBAL CLEAR INTERNAL (NON-OVERLAP MODE)- ZYLA 4.2 ONLY

Rolling Shutter Global Clear is an extended Rolling Shutter Mode. If enabled, the sensor rows are held in a global clear state to ensure there is no charge build-up on the sensor. When the camera is taken out of the Global Clear state, the exposure period begins for all rows in the sensor. When the required exposure period for Row 1 has completed, a signal frame rolling shutter readout phase begins. This means all rows have the same exposure start time but different exposure end times.

When the frame has been read out completely, the camera is returned to the Global Clear state until the user-defined cycle time is achieved or external trigger is detected (if in an external triggering mode).

Rolling Shutter Global Clear mode is best used in conjunction with illumination controlled by the Fire ALL signal.

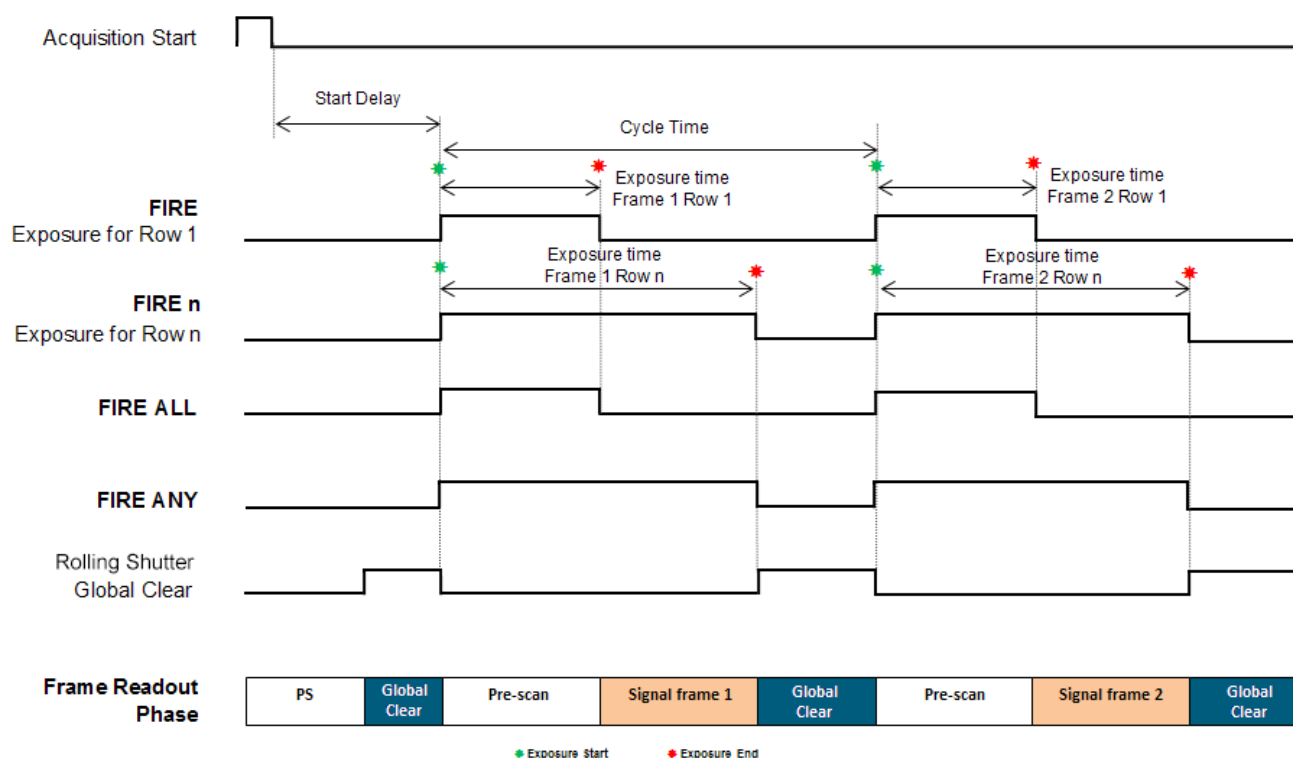


Figure 23: Rolling Shutter Global Clear- Internal (Non-Overlap Mode) - Zyla 4.2

Table 19: Rolling Shutter Global Clear Internal (Non Overlap Mode) Timing Parameters - Zyla 4.2

PARAMETER	MINIMUM	MAXIMUM
Exposure*	1 Frame + 1 Row	30s
Cycle Time (1/Frame Rate)	Exposure + 1 Frame + 5 Rows	20,000s
Acquisition Start Delay	-	2 Rows

*The exposure time reported is the exposure time for ROW 1 and the period for which FIRE ALL is HIGH.

2.6.2.10 ROLLING SHUTTER GLOBAL CLEAR EXTERNAL/SOFTWARE TRIGGERING (NON-OVERLAP MODE) - ZYLA 4.2 ONLY

In this section, both External and Software Trigger are described in the same diagram as the acquisition sequence is the same. The trigger event can either be from the EXT Trigger input or sent via software. While waiting on the trigger event, the sensor is held in a Global Clear State which ensures that charge build up on the sensor is kept to a minimum while waiting for the trigger event. The ARM signal is asserted to indicate it is ready to detect an incoming trigger input. Once the trigger event is detected, Global Clear goes LOW a signal frame read out phase begins.

When the required exposure period for Row 1 has completed, a signal frame rolling shutter readout phase begins. When the frame has been read out completely, the Global Clear is held HIGH until the next trigger event is detected.

The external trigger is fed via the EXT Trigger input on the I/O Connector on the camera head.

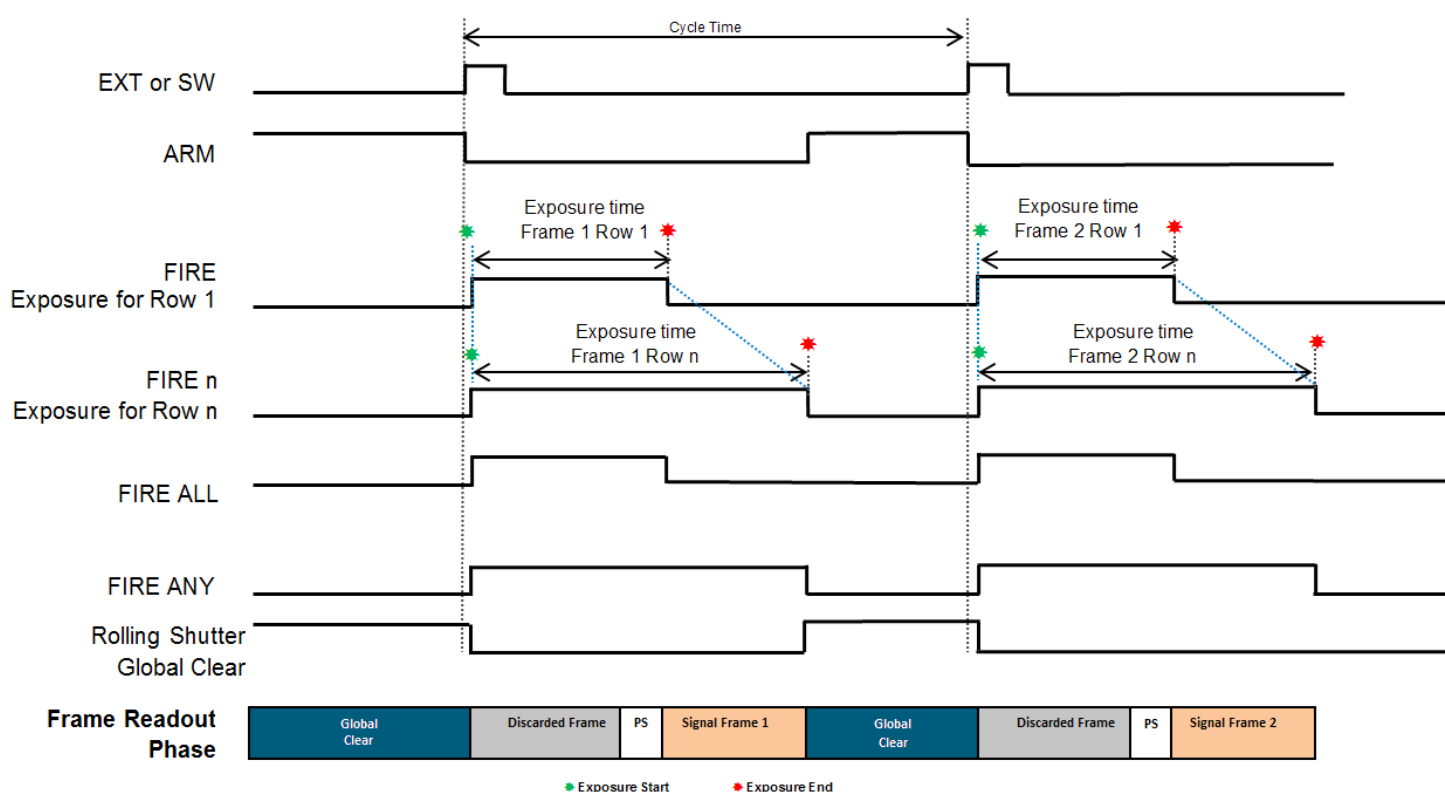


Figure 24: Rolling Shutter Global Clear External/Software Triggering (Non-Overlap Mode) - Zyla 4.2

Table 20: Rolling Shutter Global Clear External/Software Triggering (Non-Overlap Mode) Timing Parameters- Zyla 4.2

PARAMETER	MINIMUM	MAXIMUM
Exposure*	1 Frame + 1 Row	30s
Cycle Time (1/Frame Rate)	Exposure + 1 Frame + 5 Rows	-
External Start Delay	0	1 Row
EXT Trig Pulse Width	2 Sensor Speed Clock Cycles	-

*The exposure time reported is the exposure time for ROW 1 and the period for which FIRE ALL is HIGH.

2.6.2.11 ROLLING SHUTTER GLOBAL CLEAR EXTERNAL EXPOSURE TRIGGERING (NON-OVERLAP MODE) - ZYLA 4.2 ONLY

While waiting on the trigger event, the sensor is held in a Global Clear State which ensures that charge build up on the sensor is kept to a minimum while waiting for the trigger event. The ARM signal is asserted to indicate it is ready to detect an incoming trigger input.

Once the external trigger input goes HIGH, Global Clear goes LOW and the exposure begins.

When the external trigger input goes LOW, a signal frame read out phase begins. When the frame has been read out completely, the Global Clear is held HIGH until the next trigger event is detected.

The **external trigger pulse width** defines the exposure time for Row 1, and is coincident with the exposure time for Row 1. The end of exposure for Row 2 will be delayed by one row time relative to Row 1, Row 3 will be delayed by one row relative to Row 2 and so forth for the remaining rows, hence the exposure time for each row is different as all rows have the same exposure start point in time but different exposure end times.

The **period of the external trigger pulse** defines the overall cycle time. If the width of the trigger event is less than the frame read out time, the falling edge will be missed and a subsequent falling edge will be required to end the exposure.

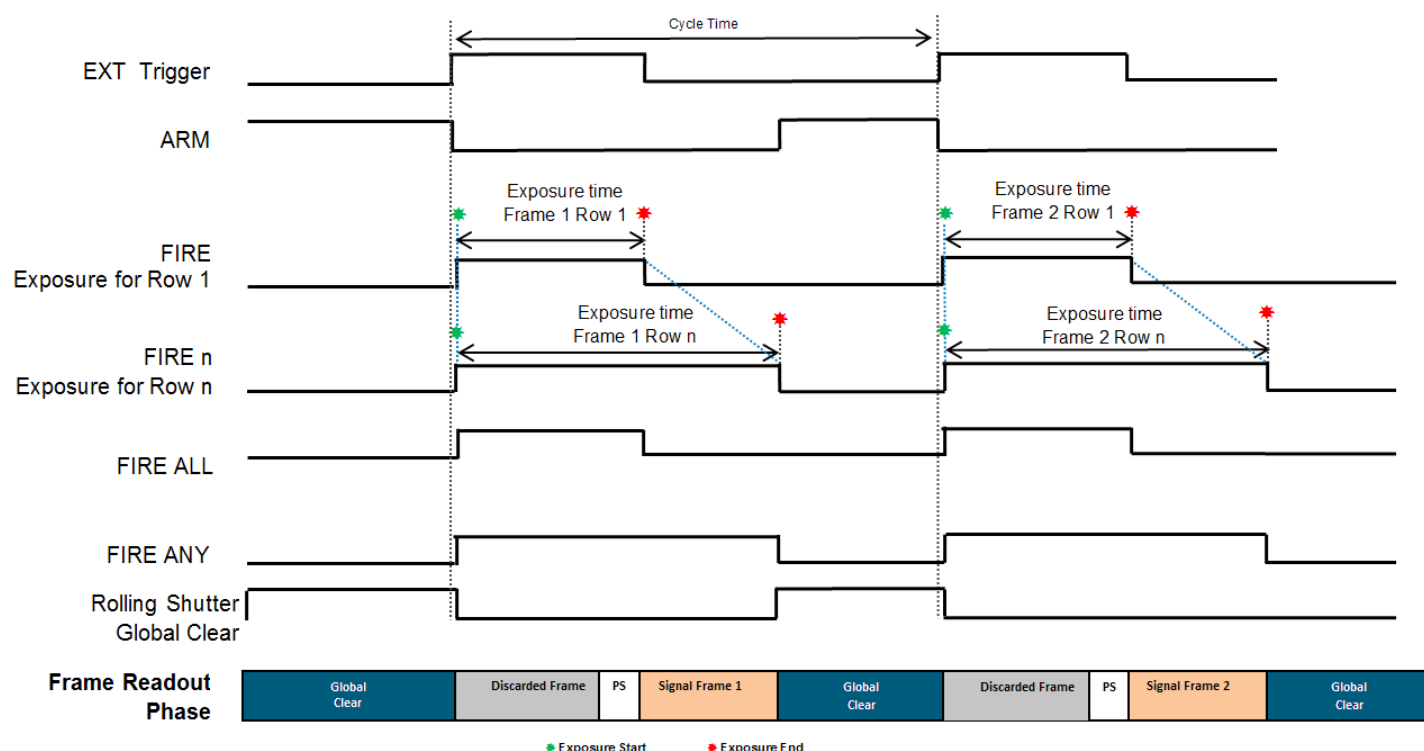


Figure 25: Rolling Shutter Global Clear External Exposure Triggering (Non-Overlap Mode) - Zyla 4.2

Table 21: Rolling Shutter External Exposure Triggering (Non Overlap Mode) Timing Parameters - Zyla 4.2

PARAMETER	MINIMUM	MAXIMUM
Exposure*	1 Frame + 1 Row	30s
Cycle Time (EXT Trig Period)	Exposure + 1 Frame + 5 Rows	-
External Start Delay	0	1 Row
EXT Trig Pulse Width	1 Frame + 1 Row	30 s

*The exposure time reported is the exposure time for ROW 1 and the period for which FIRE ALL is HIGH.

2.6.2.12 ROLLING SHUTTER TRIGGERING CONSTRAINTS

The table below shows a summary of constraints when operating in Rolling Shutter mode:

Table 22: Rolling Shutter Mode Triggering Constraints

ROLLING SHUTTER TRIGGERING MODES		EXPOSURE RANGE	MAX TRIGGER JITTER	MIN TRIGGER PULSE WIDTH	FAST EXPOSURE SWITCHING SUPPORTED
Internal (Non Overlap)	User settable exposure time. User settable Cycle Time.	1 Row to 30 s	-	-	✓
Internal Overlap (Overlap On)	User settable exposure time.	1 Row to 30 s	-	-	✓
External (Non Overlap)	Exposure time user settable. Cycle Time controlled via external trigger pulse	3 Rows to 30 s	1 Row	2 Sensor Clocks	✓
Software (Non Overlap)	Exposure time user settable. Cycle Time controlled via software trigger function	3 Rows to 30s	1 Row	2 Sensor Clocks	✓
External Exposure (Non Overlap)	Exposure Time controlled by width of external trigger pulse. Cycle Time controlled via period of external trigger pulse.	3 Rows to 30s	1 Row	2 Sensor Clocks	x
External Exposure (Overlap On)	Exposure time controlled by period of external trigger pulse	(1 Frame + 1 Row) to 30 s	1 Row	2 Sensor Clocks	x
External Start (Non Overlap)	User settable exposure time. User settable Cycle Time.	1 Row to 30 s	2 Frames 1 Row	2 Sensor Clocks	✓
Global Clear- Internal (Non Overlap)	User settable exposure time. Cycle Time is read out time for short exposures. User settable Cycle time for long exposures. Zyla 4.2 only.	(1 Frame + 1 Row) to 30 s	1 Row	-	✓
Global Clear External/ Software Triggering (Non-Overlap)	Exposure time user settable. Cycle Time controlled via period of external trigger pulse. Zyla 4.2 only.	(1 Frame + 1 Row) to 30 s	1 Row	2 Sensor Clocks	✓
Global Clear External Exposure (Non-Overlap)	Exposure Time controlled by width of external trigger pulse. Cycle Time controlled via period of external trigger pulse. Zyla 4.2 only.	(1 Frame + 1 Row) to 30 s	1 Row	1 Frame + 1 Row	✓

NOTES:

1. Exposure Time granularity for all modes is sensor row based so the exposure time in the dialog will always be rounded to the nearest integer number of row readout times.
2. 1 Row is the time taken to perform a full row readout (2624 clock cycles Zyla 5.5; 2592 clock cycles Zyla 4.2), i.e. depends on readout rate.
3. 1 Frame is the image height of a sensor half, defined in Rows and determined by the ROI configured, e.g. full frame. This is the number of rows that is read out by each sensor half. The frame readout time for any ROI can be requested via SDK function. (Refer to SDK3 User Guide for details).
4. A short exposure is referred to an exposure that is less than a Frame Readout Time. (Depends on ROI selected).
5. A long exposure is referred to an exposure time that is greater than a Frame Readout Time. (Depends on ROI selected).

2.6.3 GLOBAL SHUTTER TRIGGERING MODES (Zyla 5.5 ONLY)

Global Shutter triggering modes are available for Zyla 5.5 only. Global Shutter can also be thought of as a ‘snapshot’ exposure mode, meaning that all pixels of the array are exposed simultaneously. Before the exposure begins, all pixels in the array are cleared of charge using the Global Clear. At the start of the exposure each pixel simultaneously begins to collect charge and is allowed to do so for the duration of the exposure time. At the end of exposure, all pixels transfer their accumulated charge simultaneously to their read out node. Global Shutter requires a reference frame to be read out of the sensor in addition to the signal frame, therefore effectively halving the frame rate that would have been achieved in Rolling Shutter mode.

The Global Shutter signals shown in the diagrams are as follows:

- **Acquisition Start:** This is an internal pulse purely for illustrative purposes and indicates when the camera receives a command from software to start the pre-programmed acquisition sequence.
- **FIRE:** In Global Shutter mode, the FIRE output indicates the exposure period, which is identical for all pixels. This pulse is available to the user via the FIRE output pin.
NOTE: In Global Shutter Mode the behaviour of FIRE Row n, FIRE ALL and FIRE ANY are identical to that of FIRE and therefore not shown in the diagrams.
- **ARM:** The ARM output from the camera is used for external and software triggering modes to indicate when the camera is ready to accept another incoming trigger pulse.
- **Global Clear:** Global Shutter uses Global Clear to begin a new exposure. When this pulse is HIGH, charge is drained from every pixel thus preventing the accumulation of charge on the sensor. When the pulse is LOW, any photo-electrons generated are accumulating within the pixels, ready for transfer to the sense node for subsequent readout. The falling edge indicates the start of an exposure.
- **Charge Transfer:** This signal indicates when charge in the pixel is transferred to the measurement node and effectively ends the exposure. The charge is transferred while the pulse is HIGH and is shown in the diagrams to indicate there is a specified time between reading out the reference and signal frames.
- **Frame Read Out Phase:** This signal indicates when reference and signal frames are read out of the sensor.
- **InterFrame:** The interframe defines the minimum time taken between reference and signal frame readouts.
- * : Marks the start of an exposure.
- * : Marks the end of exposure.

The timing tables accompanying each of the triggering diagrams that follow indicate the exposure and cycle times achievable in each triggering mode. These are based on Frame and Row Periods as shown below:

Table 23: Timing Parameters based on Sensor Clock Speed for Global Shutter

PARAMETER	SENSOR READOUT RATE	
	200 MHz	560 MHz
1 Row (2624 clock cycles)	25.41 μ s	9.24 μ s
1 Full Frame (2160 rows)	27.44 ms	9.98 ms
Charge Transfer Time	5.5 μ s	2 μ s
InterFrame (9 Rows)	228.7 μ s	83.2 μ s

1 Row is the time taken to perform 2624 clock cycles. The sensor is split into 2 halves with each having an independent data output from the sensor. This means the Frame Period is 1080 rows x 2624 clock cycles.

2.6.3.1 GLOBAL SHUTTER INTERNAL TRIGGERING (NON-OVERLAP MODE)

In Internal non-overlap modes, a new exposure can not start until the previous exposure has been read out. The exact acquisition sequence depends on the exposure time. The two scenarios are shown in **Figure 26** and **Figure 27**. If the exposure time is less than the time it takes to read out a frame, then the exposure period occurs between reading out the reference and the signal frames as shown in **Figure 26**.

In this scenario, the reference frame is read out before the Global Clear is performed. The negative edge of the Global Clear begins an exposure. After the user-defined exposure time, the Charge Transfer pulse goes HIGH to transfer charge from all pixels in the sensor. The signal frame is then read out.

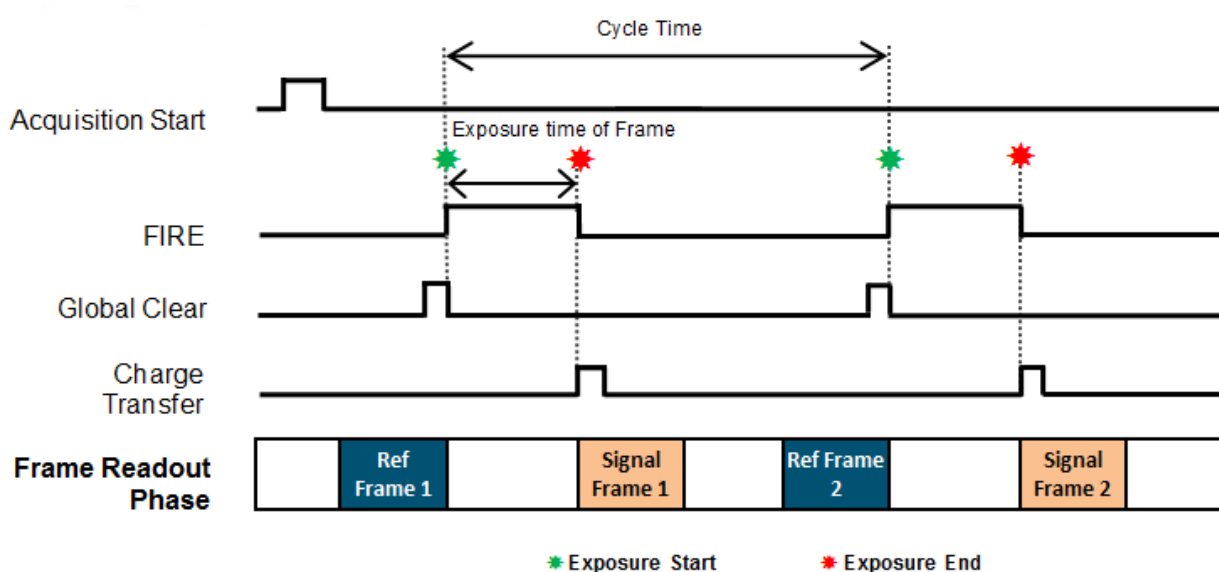


Figure 26: Global Shutter Internal Triggering – Short Exposures

Table 24: Global Shutter Internal Triggering - Short Exposures Timing Parameters

PARAMETER	MINIMUM	MAXIMUM
Exposure	1 Row	1 Frame + 3 Rows
Cycle Time (1/Frame Rate)	Exposure + 2 Frames + 2 InterFrames	Exposure + 2 Frames + 2 InterFrames
Acquisition Start Delay	0	1 Frame

If the exposure time is greater than a frame read out time, the exposure starts first by pulsing the Global Clear. The reference frame is read out during the exposure such that the end of the reference read out is coincident with the end of the exposure. The Charge Transfer pulse then goes HIGH to transfer charge from all pixels in the sensor. Finally, the signal frame is then read as shown below.

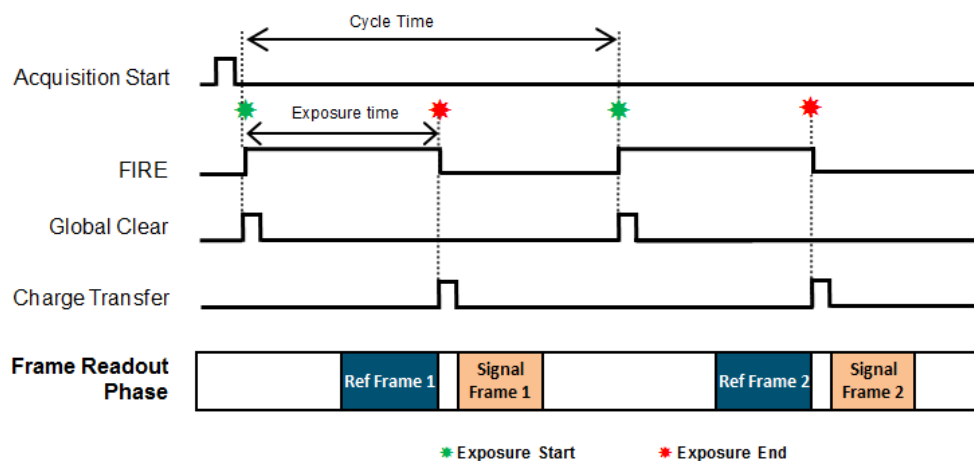


Figure 27: Global Shutter Internal Triggering – Long Exposures

Table 25: Global Shutter Internal Triggering - Long Exposures Timing Parameters

PARAMETER	MINIMUM	MAXIMUM
Exposure	1 Frame + 4 Rows	30 s
Cycle Time (1/Frame Rate)	Exposure + 1 Frames + 1 InterFrame + 6 Rows	20,000 s
Acquisition Start Delay	0	1 Row

2.6.3.2 GLOBAL SHUTTER INTERNAL TRIGGERING (OVERLAP MODE)

In Internal Triggering in Overlap Mode, the read out of an exposure overlaps with the next exposure. This allows the user to maximise the Cycle Time for a given exposure time. The absolute maximum frame rate achievable is the time taken to read out both the Reference and Signal Frame from the sensor. As in Non-Overlap Mode, the user has control over the exposure time and cycle time.

In this scenario, the exposure begins by pulsing the Global Clear. The reference frame is read out during the exposure such that the end of the reference read out is coincident with the end of the exposure. The Charge Transfer pulse then goes HIGH to transfer charge from all pixels in the sensor. The signal frame is then read out. During this read out, a new exposure begins by pulsing the Global Clear.

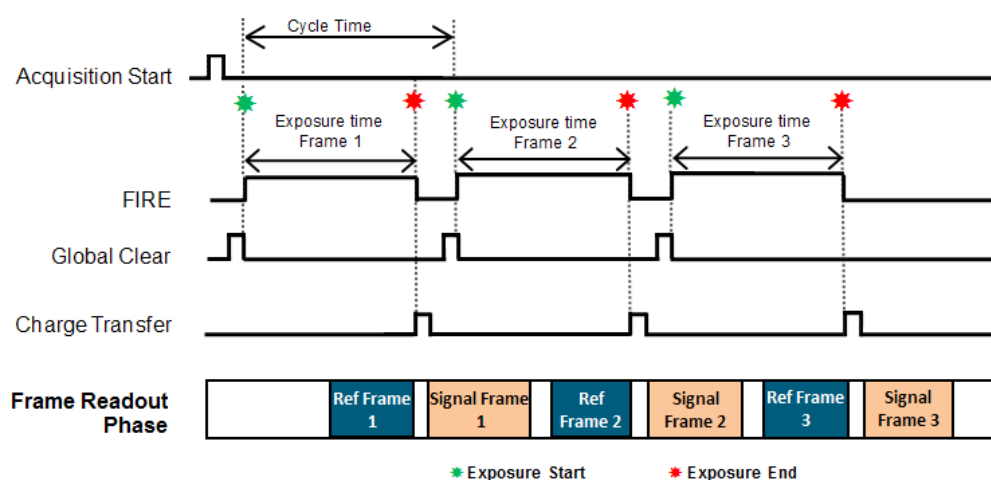


Figure 28: Global Shutter Internal Triggering (Overlap Mode)

Table 26: Global Shutter Internal Triggering (Overlap Mode) Timing Parameters - Cycle Time Dependency on Exposure

PARAMETER	MINIMUM	MAXIMUM
Exposure	1 Frame + 1 InterFrame + 1 Row	2 Frames + 1 InterFrame
Cycle Time (1/Frame Rate)	2 Frames + 2 InterFrames + 1 Row	2 Frames + 2 InterFrames + 1 Row
Acquisition Start Delay	0	1 Row
Exposure	2 Frames + 1 InterFrame + 1 Row	30 s
Cycle Time (1/Frame Rate)	Exposure + 1 InterFrame + 1 Row	30 s
Acquisition Start Delay	0	1 Row

NOTE: The table shows that the cycle time depends on the exposure selected – within the exposure range detailed in the first row the cycle time is constant, however for exposures of (2 Frames + 1 Interframe + 1 Row) or longer the cycle time increases with exposure.

2.6.3.3 GLOBAL SHUTTER EXTERNAL/SOFTWARE TRIGGERING

In this section, both External and Software Trigger are described in the same diagram as the acquisition sequence is the same. The trigger event can either from the EXT input or sent via software. While waiting on the trigger event, the sensor is put into a “pre-scan read out cycle”. On detection of the trigger, the Global Clear line is pulsed to clear the charge from the sensor. The exposure period then starts and lasts for the user defined exposure time. The reference frame is read out during the exposure such that the end of the reference read out is coincident with the end of the exposure. The Charge Transfer pulse then goes HIGH to transfer charge from all pixels in the sensor. Finally, the signal frame is then read. The camera then returns to the “pre-scan read out cycle” awaiting for the next trigger.

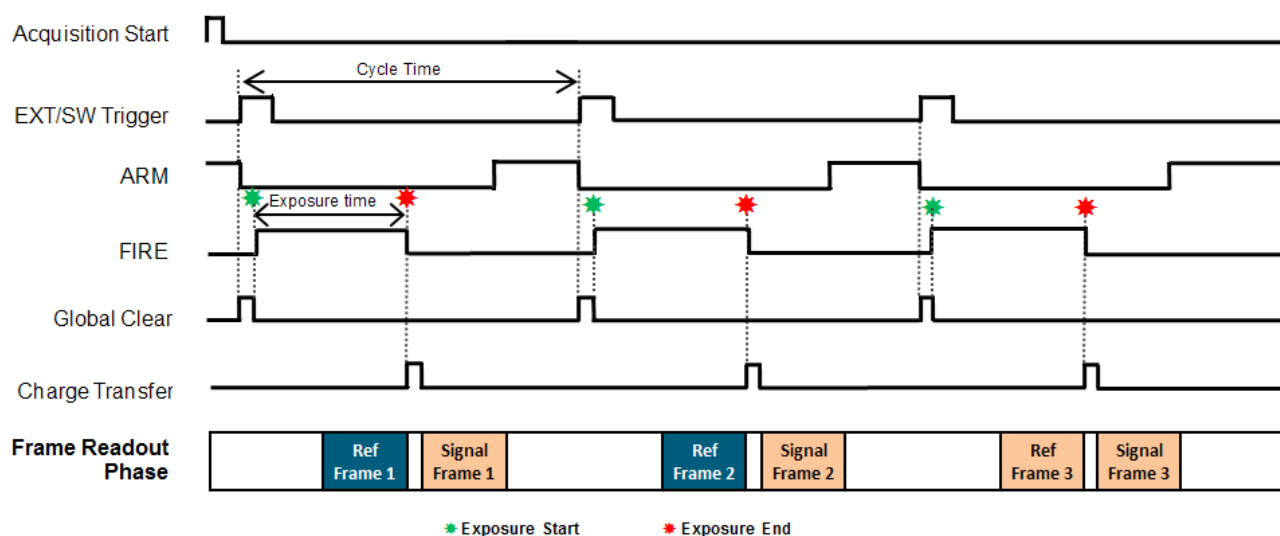


Figure 29: Global Shutter External/Software Triggering (Non Overlap Mode)

Table 27: Global Shutter External/Software Triggering Timing Parameters (Non Overlap Mode)

PARAMETER	MINIMUM	MAXIMUM
Exposure	1 Frame + 4 Rows	30 s
Cycle Time (1/Frame Rate)	Exposure + 1 Frame + 1 InterFrame + 5 Rows	
Acquisition Start Delay	1 Row	2 Rows
EXT Trig Pulse Width	2 Sensor Speed Clock Cycles	

With overlap on (**Figure 30**), the camera can accept a trigger to begin the next exposure prior to the signal frame readout completing. Note that external triggers will be missed if it is too fast- refer to **Table 28** for guidelines on time restrictions.

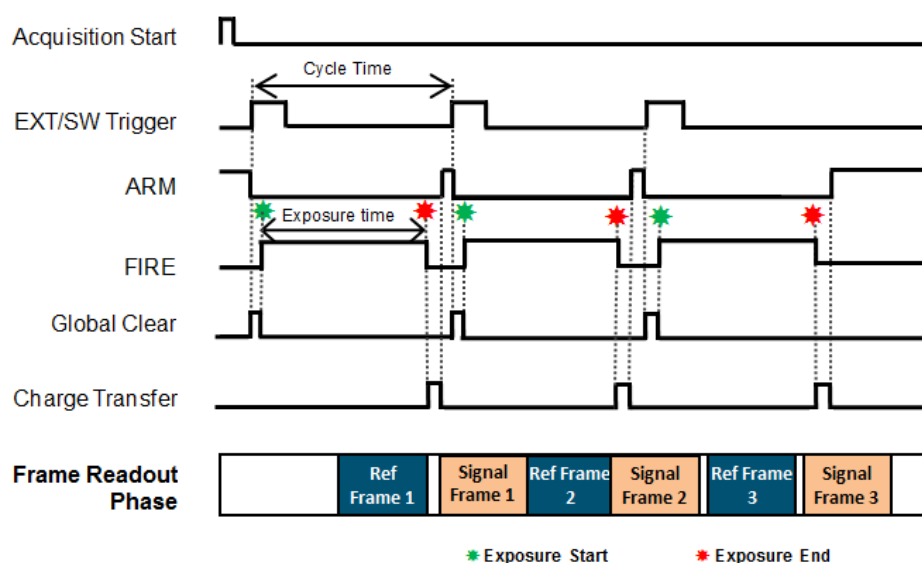


Figure 30: Global Shutter External/Software Triggering (Overlap Mode)

Table 28: Global Shutter External/Software Triggering Timing Parameters (Overlap On) - Cycle Time Dependent on Exposure

PARAMETER	MINIMUM	MAXIMUM
Exposure	1 Frame + 1 InterFrame + 1 Row	2 Frames + 1 InterFrame + 4 Rows
Cycle Time (1/Frame Rate)	2 x (1 Frame + 1 InterFrame + 1 Row)	
External Start Delay	1 Row	2 Rows
EXT Trig Pulse Width	2 Sensor Speed Clock Cycles	-
Exposure	2 Frames + 1 InterFrame + 5 Rows	30 s
Cycle Time (1/Frame Rate)	Exposure + 1 InterFrame + 2 Rows	
External Start Delay	1 Row	2 Rows
EXT Trig Pulse Width	2 Sensor Speed Clock Cycles	-

2.6.3.4 GLOBAL SHUTTER EXTERNAL EXPOSURE TRIGGERING (NON-OVERLAP MODE)

While waiting on the trigger event, the sensor is put into a “pre-scan read out cycle”. On detection of the trigger event, the Global Clear is pulsed to clear the charge from the sensor. The exposure period then starts and lasts for the width of the External Trigger. During the exposure period, a reference frame is read immediately as the system has no method to know how long the External Trigger will be HIGH for. When the external trigger input goes LOW, the exposure ends by transferring the charge in all pixels simultaneously to the measurement node. A signal frame read out then begins. The period of the external trigger pulse defines the overall cycle time. If the width of the trigger event is less than the frame read out time, the falling edge will be missed and a subsequent falling edge will be required to end the exposure.

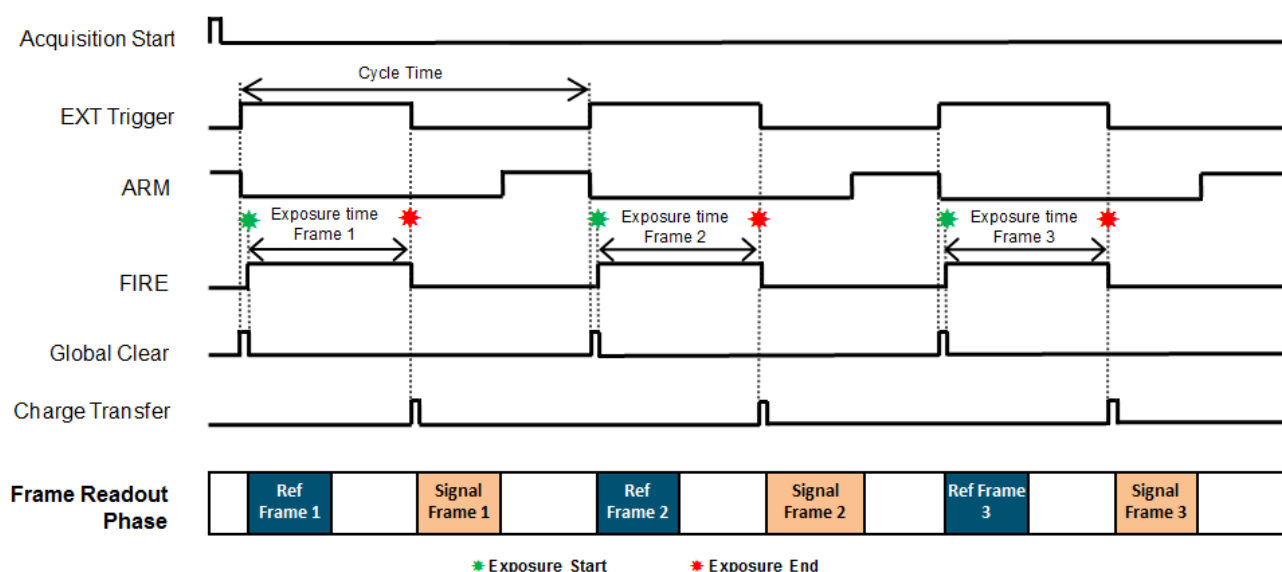


Figure 31: Global Shutter External Exposure Triggering (Non-Overlap Mode)

Table 29: Global Shutter External Exposure Triggering – Non Overlap Mode Timing Parameters

PARAMETER	MINIMUM	MAXIMUM
Exposure (EXT Trig Pulse Width)	1 Frame + 3 Rows	30 s
Cycle Time (EXT Trig Period)	EXT Trig Pulse Width + 1 Frame + 1 InterFrame + 6 Rows	
External Start Delay	1 Row	2 Rows
EXT Trig Pulse Width	1 Frame + 4 Rows	

NOTE: The table shows that the cycle time depends on the exposure selected – within the exposure range detailed in the first row the cycle time is constant, however for exposures of (2 Frames + 1 Interframe + 5 Rows) or longer the cycle time increases with exposure.

2.6.3.5 GLOBAL SHUTTER EXTERNAL EXPOSURE TRIGGERING (OVERLAP MODE)

In overlap mode, every positive edge of an External trigger will trigger a signal frame read out and start a new exposure. The period of External trigger pulse defines both the exposure time and cycle time. Note that when an acquisition starts, the first positive edge of the trigger will initiate the first exposure but also output a frame that has an incorrect exposure which is therefore discarded. The next positive edge of the trigger will end the first exposure and start a new frame read out. From the figure below, it can be seen that the minimum exposure time is approximately two frame read out periods. The first read out period is used to read out the “Signal” frame while the second is used to read out the “Reference” frame for the subsequent “Signal” frame.

If the period of the external trigger is less than the minimum period discussed above any positive edge occurring during the “Signal” and “Reference” frames will be ignored. The ARM signal indicates when positive going edges will be accepted.

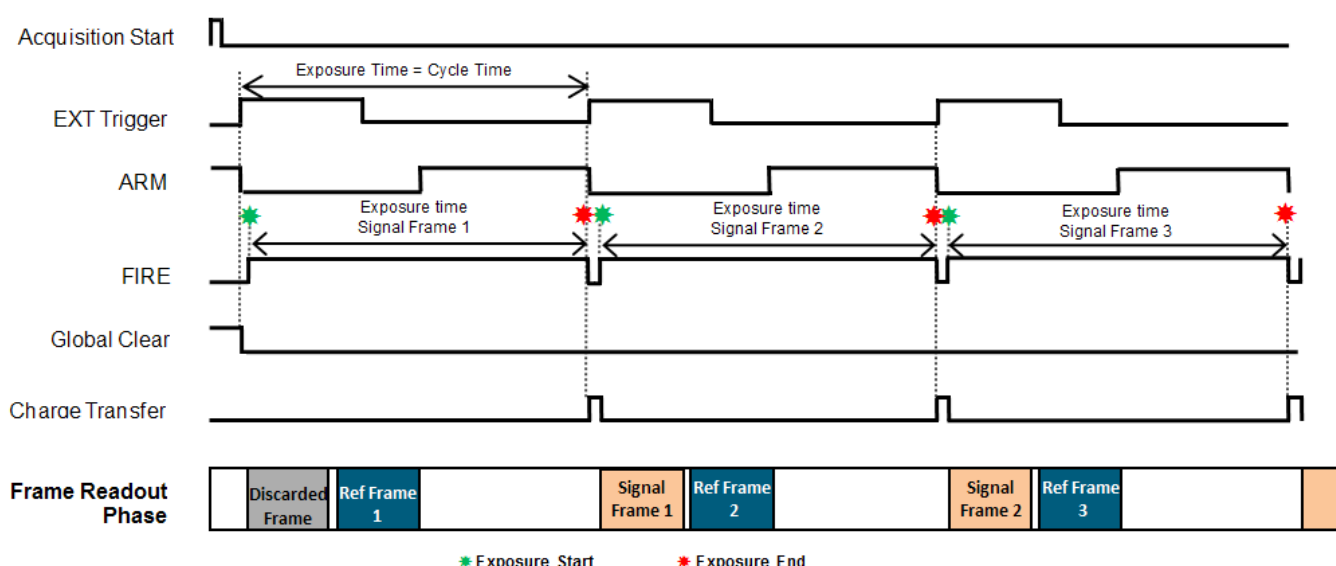


Figure 32: Global Shutter External Exposure Triggering (Overlap Mode)

Table 30: Global Shutter External Exposure (Overlap Mode) Timing Parameters

PARAMETER	MINIMUM	MAXIMUM
Exposure (EXT Trig Period)	2 Frames + 2 InterFrames	30 s
Cycle Time (EXT Trig Period)	2 Frames + 2 InterFrames	30 s
External Start Delay (1st Frame Only)	0	1 Row + Charge Transfer Time
External Start Delay (Frame #2... Frame #n)	Charge Transfer Time	Charge Transfer Time
EXT Trig Pulse Width	2 Sensor Speed Clock Cycles	-

NOTE: In global shutter external exposure mode, taking exposures in darkness of 270 ms or more will lead to increased noise. A reference frame is taken at the start of the exposure and the camera then idles until the end is signalled, which triggers the image readout. The increased noise is coming from the separation of the reference frame and the image frame. This is less prominent with deeper cooling.

2.6.3.6 GLOBAL SHUTTER EXTERNAL START TRIGGERING

In this mode the camera will wait for a single external trigger event. Once this external trigger event is detected, the camera will progress as if the camera was in internal trigger mode. The ARM signal indicates to the user when the camera is ready to detect an External Start Trigger. The diagram below shows the External Start used in an Internal Trigger Non-Overlap Mode.

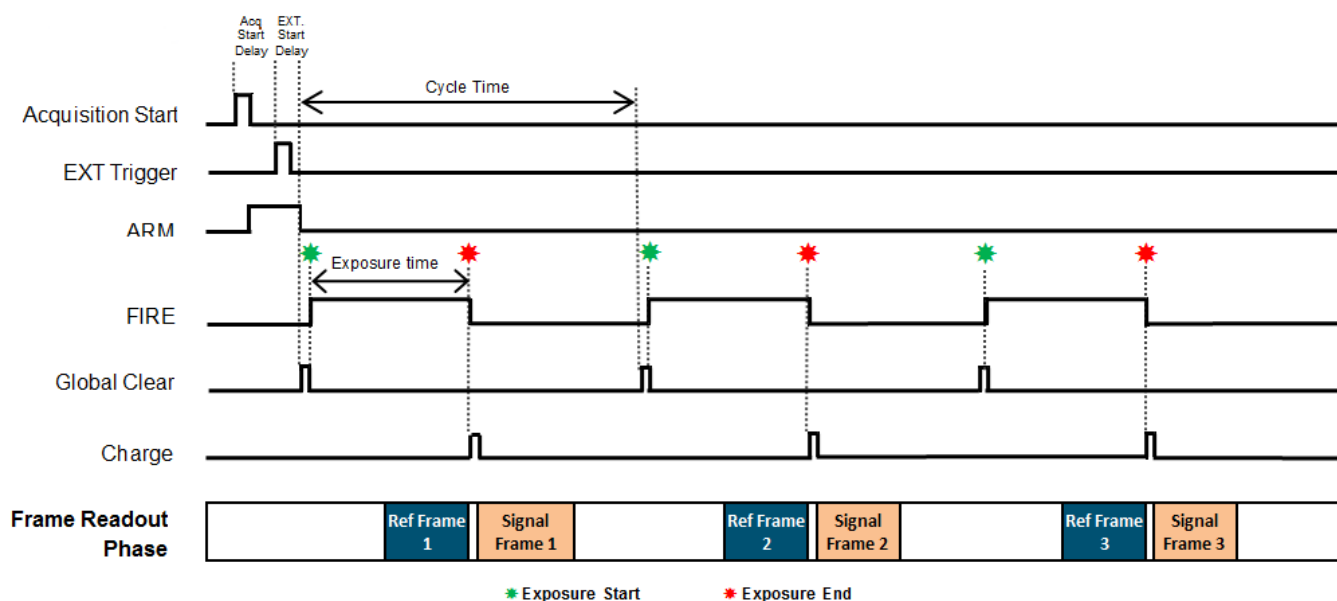


Figure 33: Global Shutter External Start Triggering

Table 31: Global Shutter External Start Triggering Timing Parameters (Overlap Off)

PARAMETER	MINIMUM	MAXIMUM
Exposure	1 Row	1 Frame + 3 Rows
External Start Delay	0	1 Frame
EXT Trig Pulse Width	2 Sensor Clock Speed Cycles	

Exposure	1 Frame + 4 Rows	30 s
External start Delay	1 Row	2 Rows
EXT Trig Pulse Width	2 Sensor Clock Speed Cycles	

Table 32: Global Shutter External Start Triggering Timing Parameters (Overlap On)

PARAMETER	MINIMUM	MAXIMUM
Exposure	1 Frame + 1 InterFrame + 1 Row	30 s
External Start Delay	1 Row	2 Rows
EXT Trigger Pulse Width	2 Sensor Clock Speed Cycles	

2.6.3.7 GLOBAL SHUTTER TRIGGERING CONSTRAINTS

The table below shows a summary of constraints when operating the Zyla 5.5 model in Global Shutter mode:

Table 33: Global Shutter Mode Triggering Constraints

GLOBAL SHUTTER TRIGGERING MODES		EXPOSURE RANGE	MAX TRIGGER JITTER	MIN TRIGGER PULSE WIDTH	FAST EXPOSURE SWITCHING SUPPORTED
Internal (Non Overlap)	User settable exposure time and Cycle Time	Short Exp: 1 Row to (1 Frame + 3 Rows)	1 Frame	-	✓
		Long Exp: (1 Frame + 4 Rows) to 30 s	1 Row	-	✓
Internal Overlap (Overlap On)	User settable exposure time	(1 Frame + 1 InterFrame + 1 Row) to 30 s	1 Row	-	✓
External/Software (Non Overlap)	User settable exposure Time. Cycle Time controlled via period of external trigger pulse	(1 Frame + 4 Rows) to 30 s	1 Row	2 Sensor Clocks	✓
External/Software (Overlap On)	User settable exposure time. Cycle Time controlled via software trigger function	(1 Frame + 1 InterFrame + 1 Row) to 30 s	1 Row	2 Sensor Clocks	✓
External Exposure (Non Overlap)	Exposure Time controlled by width of external trigger pulse. Cycle Time controlled via period of external trigger pulse.	(1 Frame + 3 Rows) to 30 s	1 Row	1 Frame + 3 Rows	✗
External Exposure (Overlap On)	Exposure time controlled by period of +ve edge of external trigger pulses. Cycle Time is exposure time	(2 Frames + 2 InterFrames) to 30 s	1 Row	2 Sensor Clocks	✗
External Start (Non Overlap)	User settable exposure time and Cycle Time but sequence initialized via external trigger pulse	Short Exp: 1 Row to (1 Frame + 3 Rows) Long Exp: (1 Frame + 4 Rows) to 30 s	1 Frame 1 Row	2 Sensor Clocks	✓ ✓
External Start (Overlap On)	User settable exposure time	(1 Frame + 1 InterFrame + 1 Row) + 30 s	1 Row	2 Sensor Clocks	✓

NOTES:

1. Exposure Time granularity for all modes is sensor row based so the exposure time in the dialog will always be rounded to the nearest integer number of row readout times.
2. 1 Row is the time taken to perform a full row readout (2624 clock cycles), i.e. depends on readout rate.
3. 1 Frame is the image height of a sensor half, defined in Rows and determined by the ROI configured, e.g. full frame = 1080. This is the number of rows that is read out by each sensor half. The frame readout time for any ROI can be requested via SDK function. (Refer to SDK3 User Guide for details).
4. A short exposure is referred to an exposure that is less than a Frame Readout Time. (Depends on ROI selected).
5. A long exposure is referred to an exposure time that is greater than a Frame Readout Time. (Depends on ROI selected).
6. Image acquisition must be stopped to change between short and long exposures. Fast Exposure Switching can occur in the triggering modes shown above but the exposure time is then limited to the exposure range shown for each sequence.

2.7 ACQUISITION MODES

The following acquisition modes can be supported:

- Single Scan
- Kinetic Series
- Accumulate
- Run Till Abort

NOTES:

1. The term 'User Frame', in this section refers to a single frame in Rolling Shutter mode and a reference/image frame pair in Global Shutter mode.
2. The term 'valid trigger' refers to a trigger that is applied when the camera is ready to accept it.

2.7.1 SINGLE SCAN

Single Scan refers to an acquisition in which only one user frame is transmitted from the camera.

A user frame is output from the sensor on receipt of a valid trigger of the selected type and then transmitted from the camera. Note that any subsequent triggers within the same acquisition are ignored.

2.7.2 KINETIC SERIES

Kinetic Series refers to an acquisition in which a finite number of user frames are transmitted from the camera. The number of frames in a Kinetic Series is defined by the user. One user frame is output from the sensor on receipt of each valid trigger of the selected type. Valid triggers continue to output user frames from the sensor until the defined number of user frames has been reached. Frames are transmitted from the camera, depending upon the connectivity i.e. 3-tap or 10-tap Camera Link. Note that after the required number of valid triggers has been received, any subsequent triggers within the same acquisition are ignored.

2.7.3 ACCUMULATE

Accumulate refers to an acquisition in which a number of frames in a series are accumulated together into a single image. This accumulation of user frames is performed off-camera. Either all the user frames in a series are accumulated to give a single accumulated image or a smaller number of user frames in the series are accumulated to give a series of accumulated images.

2.7.4 RUN TILL ABORT ACQUISITION

Run Till Abort refers to an acquisition in which an infinite number of user frames can be transmitted from the camera and the acquisition will continue to run until it is aborted.

One user frame will be output from the sensor on receipt of each valid trigger of the selected type. All valid triggers will output another user frame from the sensor. Frames are transmitted from the camera in as quick succession as possible.

2.7.4.1 LIVE MODE

Live Mode refers to a version of **Run Till Abort** in which each user frame will be the latest frame output by the sensor and will have the minimum amount of latency through the camera as possible.

Live mode requires the use of SW Trigger. In order to ensure that frames are buffered in the on-camera memory for the shortest possible time, the next SW trigger must not be sent until the user frame has been transmitted from the camera. This ensures that the on-camera memory only stores a single user frame at a time and no additional latency builds up.

NOTE: The frame rate achievable in Live Mode is dependent on the performance of the system that the camera is attached to.

2.7.5 FAST EXPOSURE SWITCH

During an acquisition the user can change the exposure time, within allowable limits. Once a new exposure value has been written, it will be applied to the next user frame after the current frame exposure has completed. The exposure time can be changed any number of times before the acquisition finishes

2.7.6 FRAME RATE CONTROL

If Internal Trigger is being used, the camera will trigger the sensor at the fastest possible rate by default. The user can reduce this trigger rate by defining a Frame Rate that is less than the maximum possible rate. Frame Rate must be set before the acquisition starts.

Note: This is currently only applicable to 'long exposures'.

SECTION 3: INSTALLATION

3.1 SAFETY CONSIDERATIONS

- Prior to commencing installation, please refer to the Specifications in **Section 1.4** to ensure all requirements have been met
- As part of the safety features of the product the Zyla must be powered from an SELV (safety extra low voltage) source. In the case of the specified power supply (SW3605) this unit provides an SELV output and is designed to have a protective earth connected via the earth pin on the mains plug of the unit. It is important to ensure that this is connected to the buildings protective earth system
- The equipment should be positioned so that the mains supply plug/cord can be easily accessed for disconnection

3.2 CAMERA MOUNTING

To prevent damage to the Zyla, other equipment and/or personnel the camera must be securely mounted before use. This can be using either one of the following options:

- The C-Mount, CS-Mount or F-Mount on the front face of the camera (please refer to **Section 4.2**)
- The ¼ -20 UNC threaded holes located on each side of the unit (please refer to **Appendix A**).
- The 4-off M4 threaded holes on the front face of the camera (please refer to **Appendix A**).

How to use the lens mount adaptors is shown in **Section 4.2**.

3.3 COOLANT HOSE CONNECTION AND DISCONNECTION (WATER COOLED MODELS)



CAUTION:

- Before attempting to insert or remove the coolant hose connections, ensure that all coolant has been drained from the hoses and integral coolant channel within the camera head.
- Care must be taken to avoid permanent damage to the camera system resulting from either leakage of coolant during connection / removal of hoses or spillage of any residual coolant contained within the camera head once the hoses have been removed.
- Always ensure that the temperature of the liquid coolant circulated through the camera head is above the dew point of the camera ambient temperature and humidity conditions. Refer to the Dew Point graph in **Appendix B** for guidance.
- Use of coolant at or below the dew point can result in permanent damage to the camera head, due to formation of condensation on internal components.
- Never use damaged, split or worn hoses.

Coolant Hose Connectors: Two barbed coolant hose inserts are supplied as standard with the water cooled Zyla models, suitable for connection to 6 mm (0.25") internal diameter soft PVC tubing / hose.

Recommended tubing: 10 mm (0.4") outside diameter, i.e. a wall thickness of 2 mm (0.08"). Alternative hose dimensions and materials should be thoroughly tested to ensure a leak tight seal is achieved with the barbed inserts.

3.3.1 COOLANT RECOMMENDATIONS

It is recommended that de-ionized water (without additives) is used as the coolant to prevent deposits forming. Some mains supply water is heavily mineralized (i.e. "Hard") which could cause deposits in the water circuit inside the camera. This can reduce the flow-rate and cooling efficiency.

The specified cooling performance of the camera can be achieved with coolant flow rates of 2 litres per minute, the maximum recommended pressure of coolant circulating through the camera head is 2 bar (30 PSI).

In the event that replacement hose inserts / barbs are required, please contact your local Andor representative.

3.3.2 CONNECTING THE COOLANT HOSES

1. Press the hose connector into the connection on the camera head, ensure it clicks into place and repeat for the second hose.
2. Confirm the hoses are connected securely by applying pressure on the top front of the camera body and pulling backwards on the hoses.
3. The coolant hoses are now connected.

3.3.3 DISCONNECTING THE COOLANT HOSES

1. Press the latch on the camera hose connection away from the hose.
2. Hold the latch in and pull the hose backwards.
3. The hose should release from the camera connection with little resistance.

NOTE: If the hose does not release, ensure that the latch on the camera connection is pressed in fully.

3.4 RECOMMENDED COMPUTER REQUIREMENTS

- 2.68 GHz Quad Core + 4 GB RAM
- Hard Drive:
 - Minimum 250 MB/s continuous write for 'Zyla 5.5' 3-tap model
 - Minimum 850 MB/s continuous write for 'Zyla 5.5 and Zyla 4.2' 10-tap models
 - Minimum 450 MB/s continuous write for Zyla USB 3.0 model
- PCI Express x4 or greater for 'Zyla 5.5' 3-tap
- PCI Express x8 or greater for 'Zyla 5.5 and Zyla 4.2' 10-tap model
- PCI Express x4 or greater for Zyla USB 3.0 models
- Windows (Vista, 7 and 8) or Linux

3.5 INSTALLING SOFTWARE (SOLIS OR SDK3)

NOTE: You must have administrator access on your PC to perform this installation.

The same instructions cover the installation procedure for Andor's Solis software or Andor's SDK, which is used in conjunction with third party software. If you are planning to run your camera through a third party interface you will require the Andor Drivers, called SDK3.

Switch on the PC, insert CD and run the "setup.exe" file on the CD or download location.

- Confirm the version of software.
- Follow the on screen prompts.
- Select the installation directory when prompted.
- When prompted for camera type select Zyla sCMOS.
- On the final window click on "Install". During the installation a number of other windows will appear as various drivers are installed.
- Click on the finish button when prompted.






Figure 34: Selecting the Camera Type

For more information visit: andor.com/software/software_support/

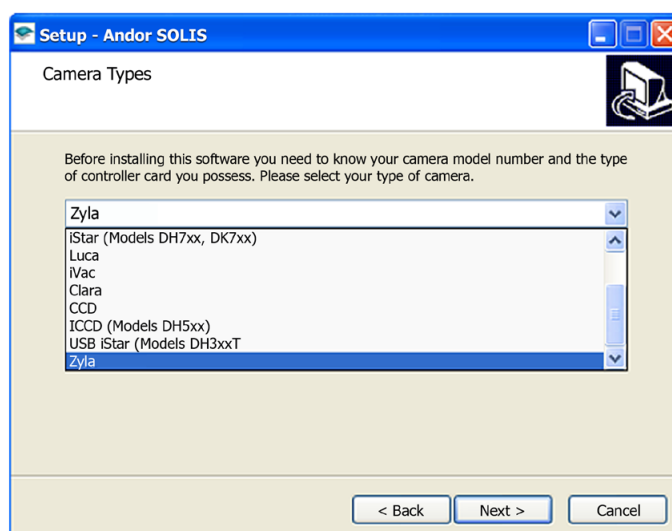3.6 INSTALLING THE CAMERA FRAMEGRABBER CARD




Ensure your PC is powered off. Ensure anti-static precautions using anti-static strap provided. Keep all screws safe.

1. Unplug all cables from the rear of the computer.
2. Remove the PC enclosure so as to gain access to the expansion slots.
3. Locate a PCIe x16, x8 or x4 slot.
 - **10-tap version:** Karbon Framegrabber card operates in x8 or greater (i.e. x16 and x8 slots) ONLY.
 - **3-tap version:** Neon Framegrabber card operates in x4 or greater (i.e. x16, x8 or x4 slots).
 - **USB 3.0 version:** Framegrabber card operates in x4 or greater (i.e. x16, x8 or x4 slots).

If you are unsure which PC slot is correct, please consult the PC user manual.

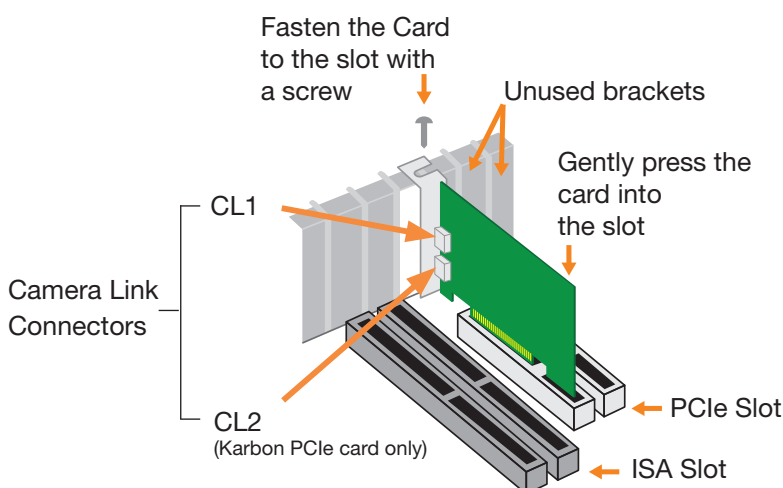4. Remove the filler bracket corresponding to the slot you intend to use.
5. Remove the controller card carefully from its protective ESD packaging and firmly press the card connector into the chosen expansion slot and ensure it is securely locked in.
6. Making sure that the card's mounting bracket is flush with any other mounting brackets or filler brackets to either side of it, secure the controller card in place.
7. Replace the cover of the computer and secure it with the mounting screws if applicable.
8. Reconnect any accessories you were using previously.

Camera Link Models



USB 3.0 Models

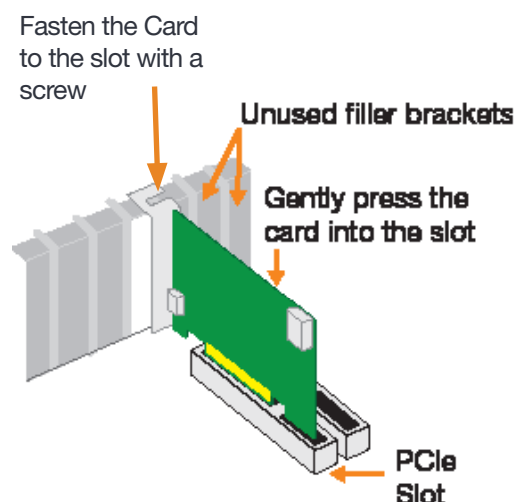


Figure 35: Installing the FrameGrabber PCIe card (Left) Camera Link versions and (Right) for USB 3.0.

3.7 INSTALLING THE PCIe CARD DEVICE DRIVER (CAMERA LINK MODELS)



Restart your PC



If you are running Windows 8, the drivers will be installed automatically during start-up. Continue to the next section.



If you are running Windows 7, the drivers will be installed automatically during start-up. Continue to the next section.



If you are running Windows Vista you may be asked to specify the directory containing the Bitflow drivers. These are located in 'C:\Bitflow SDK 5.60\PlugAndPlay'.



If you are running Windows XP, the 'Found New Hardware Wizard' dialog will be displayed following boot up. When prompted 'Can Windows connect to Windows Update to search for software?', select 'No' and click on Next. Select 'Install the software automatically' option and click on 'Install'. The Camera Link drivers will now be installed.

3.7.1 IMPORTANT INSTALLATION INFORMATION FOR USB 3.0 CAMERAS

- For USB 3.0 cameras, the drivers for the PCIe card must be installed manually by the user as they are not part of the SDK installation. Refer to the instructions on the driver CD supplied with the USB 3.0 Card.
- Power management in the Renesas USB 3.0 Host Controller must be disabled. To do this:
 - Navigate to 'Renesas USB 3.0 Host Controller' properties menu in **Device Manager** (Control Panel>Device Manager> Universal Serial Bus Controllers).
 - Select the **Advanced** tab.
 - Check the **Disable USB 3.0 power management functions** checkbox as shown in Figure 36.

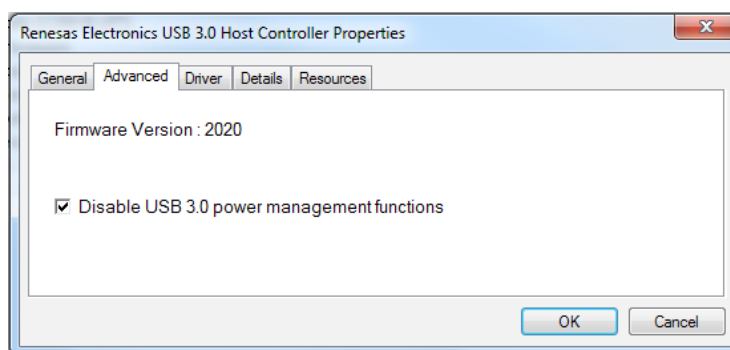


Figure 36: Disabling the power management function of the USB 3.0 Host Controller

3.8 CHECKING & SETTING BIOS OPTIONS (FOR PCs NOT SUPPLIED BY ANDOR)

Enter the BIOS menu when starting PC. For Dell workstations, press F12 at start-up and select System Setup in the One Time Boot Menu. For Dell workstations 3 options in the Performance menu of the BIOS need to be checked/set:

- C-States Control – Disable C-States
- Intel Speed-step – Disable Speed-step
- Memory Node Interleaving – Set from NUMA to SMP. Note: This option is only available on larger workstations with 2 physical processors and may have a different name- ensure that NUMA is disabled.

SECTION 4: OPERATION

4.1 CONNECTING YOUR CAMERA AND SWITCHING IT ON



Ensure that the Camera on / off switch is in the **OFF** position **AND** your PC is switched **OFF**.



For water cooled models, ensure the cooling water supply is connected prior to powering the camera (refer to Section 3.3).

- Before inserting the 12V DC power connector to the rear of the camera, ensure the orientation is correct. **Never** forcibly insert the connector. For water-cooled models, connect water cooling hoses (see Section 3.3.2) and ensure re-circulator/chiller is on.
- Securely connect the Camera to PC cable between the appropriate connector(s) on the rear of the camera and the corresponding port on the PCIe card.
(See Figure 37, please refer also to diagram in Section 3.5 for correct PCIe slot connection).

Notes:



Andor Camera Link cables are labelled to assist correct installation.



For USB 3.0 models, use only the supplied USB 3.0 cable to ensure performance.

- If required connect the I/O expander cable to the 15 way I/O connector.
- Switch the camera on using the ON/OFF switch. You should hear an audible confirmation (a short beep followed by a long beep)
- Turn on the PC.
- Start your software.
- Start acquiring data.

Camera Link Model



USB 3.0 Model



1. I/O 15-Way D Type Connector
2. CL1 connector
3. CL2 connector (10-tap cameralink versions only)
4. USB 3.0 Connector

Figure 37: Connect the Zyla to the port on the PC by the connections shown above (Left) Camera Link model, (Right) USB 3.0 model

4.2 USING THE LENS MOUNT ADAPTORS

The Zyla has a T-mount thread as standard. There are 3 standard lens mount adaptors available that can be specified at the time of ordering, enabling the Zyla to suit standard mounting options. These are as follows:

- C-mount
- CS-mount
- F-mount

The Zyla is supplied with the selected lens mount adaptor locked into a **nominal position of 1 mm** from the top surface of the Zyla faceplate (see **Figure 38**). Adaptors can be removed, inserted and adjusted (for focus optimization) as required.

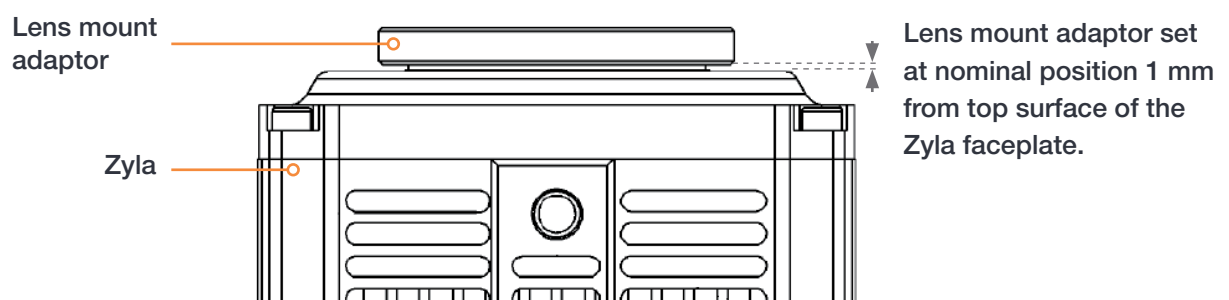


Figure 38: Lens mount adaptor fitted to Zyla (In this example a CS-mount adaptor is shown).

4.2.1 C-MOUNT ADAPTOR

The C-mount adaptor can be adjusted as follows:

1. Loosen the 2 locking screws located in the front face of the adaptor if required using supplied 1.5 mm Allen key.

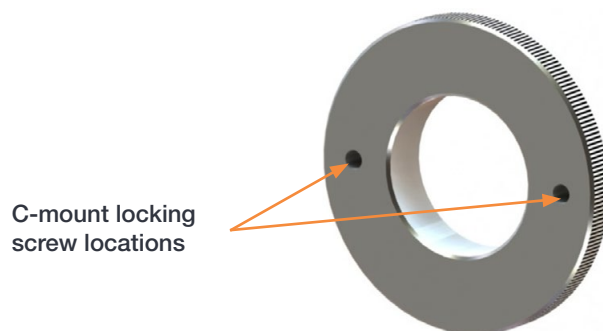


Figure 39: C-mount adaptor

2. Manually screw the C-mount into position until the desired focus is achieved. It must not be screwed the entire way into the faceplate.
3. Lock the C-mount into position using the 2 locking screws and 1.5 mm Allen key. **NOTE: Maximum torque applied must not exceed: 0.6 Nm**

4.2.2 CS-MOUNT ADAPTOR

The CS-mount adaptor can be adjusted as follows:

1. Loosen the 2 locking screws located in the front face of the adaptor if required using supplied 1.5 mm Allen key.

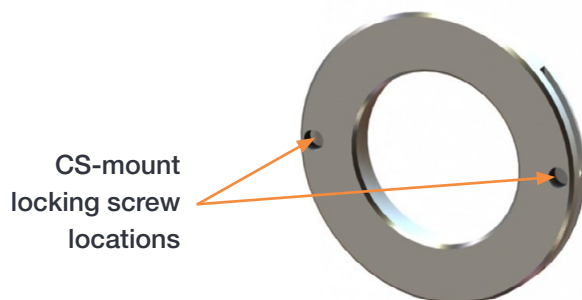


Figure 40: CS-mount adaptor

2. Manually screw the CS-mount into position until the desired focus is achieved. It must not be screwed the entire way into the faceplate.
3. Lock the CS-mount into position using the 2 locking screws and 1.5 mm Allen key. **NOTE: Maximum torque applied must not exceed: 0.6 Nm**

4.2.3 F-MOUNT ADAPTOR

The F-mount should be screwed into the camera faceplate by hand.



Figure 41: F-mount adaptor

1. Manually screw the F-mount adaptor until the desired focal position is achieved.
2. Manually tighten the collar down to the camera faceplate to lock the mount in position.

4.3 COOLING

4.3.1 INTEGRAL THERMAL PROTECTION

Care should be taken to ensure that the camera does not overheat, as this can cause system failure. Overheating may occur if either of the following situations arises:

Air-cooled Models

- The air vents on the sides of the head are accidentally blocked
- The ambient air temperature is more than 30°C (Zyla 5.5*), or 27°C (Zyla 4.2)

*For Zyla 5.5 models manufactured prior to February 2015 max. ambient is 35°C.

To protect the camera from overheating, a thermal switch has been attached to the heat sink. If the temperature of the heat sink rises above predefined limit then the current supply to the cooler will cut out and a buzzer will sound. To reset this, it is necessary to exit the software and switch off the camera, correct the issue, let the system cool then restart the camera and software.

Water-cooled Models

- The coolant system is not operating correctly

To protect the camera from overheating, a thermal switch has been attached to the heat sink. If the temperature of the heat rises above predefined limit then a buzzer will sound. To reset this, it is necessary to exit the software and switch off the camera, correct the issue, let the system cool then restart the camera and software.

4.3.2 TE COOLING (AIR-COOLED MODELS)

The Zyla sCMOS sensor is cooled using a thermoelectric (TE) cooler. TE coolers are small, electrically powered devices with no moving parts, making them reliable and convenient. A TE cooler is actually a heat pump, i.e. it achieves a temperature difference by transferring heat from its 'cold side' (the CMOS-chip) to its 'hot side' (the built-in heat sink). Therefore the minimum absolute operating temperature of the Zyla sensor depends on the temperature of the heat sink. The maximum temperature difference that a TE device can attain is dependent on the following factors:

- Heat load created by the Zyla sensor (dependant on pixel readout speed)
- Number of cooling stages of the TE cooler
- Operating current
- Hot side temperature of thermoelectric cooler

Even with a fan, a heat sink typically needs to be at least 10°C hotter than the air (room) temperature to transfer heat efficiently to the surrounding air. Therefore during normal operation the temperature of the heatsink within the camera head will be warmer than the ambient (room) temperature.

4.3.3 TE COOLING (WATER COOLED MODELS)

For water cooled models the dissipation of heat from the TE cooler is achieved by circulating coolant liquid through the heat sink. The high specific heat capacity of water provides efficient heat dissipation, allowing the sensor to be operated at -10°C. Benefits of water cooling include:

1. Reduced vibration as there is no cooling fan required, which may be useful for ultra-sensitive measurements and
2. Improved dark current.

4.4 SOFTWARE SUPPORT

The following software applications include support for the Zyla camera, check with your software supplier for applications not listed below:

- Andor Solis
- Andor SDK3
- Andor iQ
- MetaMorph
- NIS-Elements
- MicroManager

Please refer to the application Help for details on how to control any Zyla specific functionality. Andor SDK3 documentation is provided in the **Andor Software Development Kit 3.pdf** file that comes with the installation package if this has been ordered as an option.

SECTION 5: MAINTENANCE



THERE ARE NO USER-SERVICEABLE PARTS INSIDE THE CAMERA. DAMAGE CAUSED BY UNAUTHORISED MAINTENANCE OR PROCEDURES WILL INVALIDATE THE WARRANTY.

5.1 REGULAR CHECKS

- The state of the product should be checked regularly, especially the integrity of the External Power Supply and the mains cable.
- Do not use equipment that is damaged

5.2 ANNUAL ELECTRICAL SAFETY CHECKS

- It is advisable to check the integrity of the insulation and protective earth of the AC/DC converter on an annual basis, e.g. U.K. PAT testing
- Do not use equipment that is damaged

5.3 FUSE REPLACEMENT

In the U.K, Ireland and some other countries, the supplied mains cable has a BS 1363 (or Type G) plug that includes an integrated fuse. Only replace with fuse of the same type and rating for continued protection. The characteristics of a replacement fuse are as follows:

- **Rated Current:** 5 A
- **Rated Voltage:** 240 VAC
- **Size:** ¼ × 1" (6.3 × 25.4 mm) cartridge
- **Type:** BS 1362

5.4 COOLING HOSES AND CONNECTIONS

The user should routinely check all coolant hoses and connections for signs of leakage, damage or wear. All seals must be intact before powering on camera system and any worn / damaged items must be replaced immediately.

SECTION 6: TROUBLESHOOTING

This section provides useful information and solutions for some troubleshooting scenarios. If you have an issue that you are unable to rectify using this section, please contact Andor Technical Support for further advice.

6.1 CAMERA BUZZER DOES NOT SOUND ON START-UP

1. The camera buzzer should be audible momentarily when the camera is switched on.
2. If this does not occur, ensure that power is connected to the camera and the On/Off switch is set to On

6.2 CAMERA IS NOT RECOGNIZED BY PC

1. Ensure the camera is switched on
2. Check the Camera to PC connection. For Camera Link models, check that the cable is connected between the **CL1** connector position and the PCIe card in the PC. (For 10-tap variants, check that the **CL2** cable is connected). For USB 3.0 models, check that USB 3.0 cable is connected between the camera and the USB 3.0 slot on the PC.
3. Check that the frame grabber card and drivers have been correctly installed (see **Sections 3.4 and 3.5**)

6.3 BUZZER SOUNDS CONTINUOUSLY

This indicates that an “over temperature” condition has occurred within the camera. Follow the instructions below to rectify this situation:

1. Power the camera off and allow it to cool down
2. For air-cooled models, ensure fan vents are not obstructed
For water-cooled models, ensure that water coolant supply is connected and that chiller/re-circulator is operational
3. Check the camera is operating within the specified environmental conditions (see **Section 1.4**)

6.4 FAN NOT OPERATING AS EXPECTED

- To protect the internal electronics, the fan defaults to full speed if the camera heat-sink temperature exceeds 50°C.

NOTE: The fan will continue to run even if the user has switched it off via software, until the correct heat-sink temperature has been reached.

6.5 CAMERA DOES NOT COOL TO THE REQUIRED TEMPERATURE

1. Check that the operating ambient temperature is within allowable limits (see **Table 1**) when cooling the sensor to 0°C.
2. Check that the camera vents are not blocked and have sufficient clearance to allow air flow.
3. Check that the fan is switched on (or for water cooled modesl, the coolant system is opertaing correctly).

6.6 FIRE, AUX_OUT AND ARM OUTPUTS NOT FUNCTIONING CORRECTLY

These are 5V TTL outputs which should not be used to drive low impedance loads. Please refer to **Section 1.7.2** and **1.7.3** for suggested terminations.

6.7 EXTERNAL TRIGGER INPUT NOT FUNCTIONING CORRECTLY

This is a 5V TTL input which should be driven from a 5V TTL compatible source. Please refer to **Section 1.7.2** and **1.7.3** for further details.

6.8 PREVENTING CONDENSATION

NEVER USE WATER THAT HAS BEEN CHILLED BELOW THE DEW POINT OF THE AMBIENT ENVIRONMENT TO COOL THE CAMERA.

You may see condensation on the outside of the camera body if the cooling water is at too low a temperature or if the water flow is too high. The first signs of condensation will usually be visible around the connectors where the water tubes are attached. If this occurs carry out the following actions:

1. Switch off the system
2. Wipe the camera with a soft, dry cloth.

NOTE: It is likely there will already be condensation on the cooling block and cooling fins inside the camera.

3. Set the camera aside to dry for several hours before you attempt reuse.
4. Before reuse blow dry gas through the cooling slits on the side of the camera to remove any residual moisture.

Use warmer water or reduce the flow of water when you start using the device again.

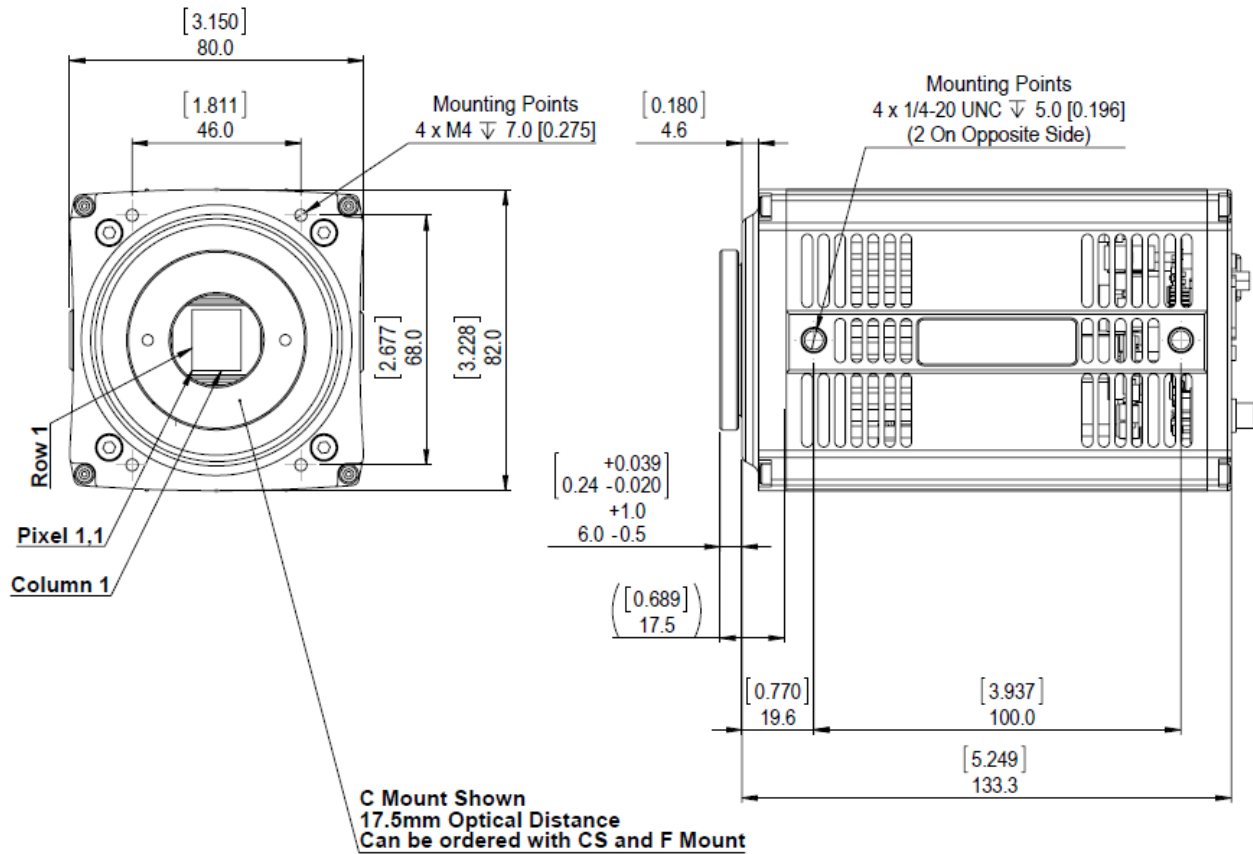
NOTE: This is not an issue when using a Re-circulator which eliminates the dew point problem.

Refer to **Appendix B** for a Dew Point Graph.

APPENDIX A: MECHANICAL DRAWINGS

Zyla Mechanical Drawings


Air Cooled Models



Zyla 5.5 3-tap version



Zyla 5.5 and Zyla 4.2 10-tap

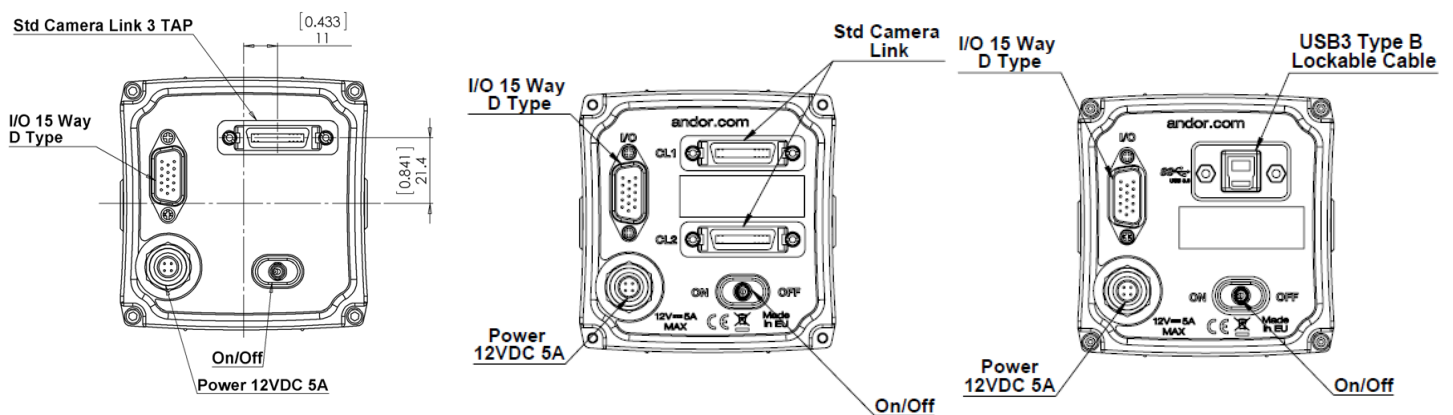Zyla 5.5 and 4.2 USB 3



Appendix

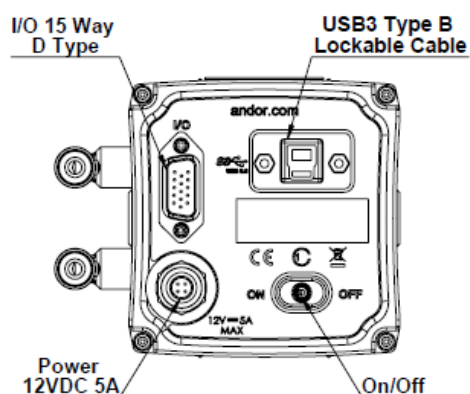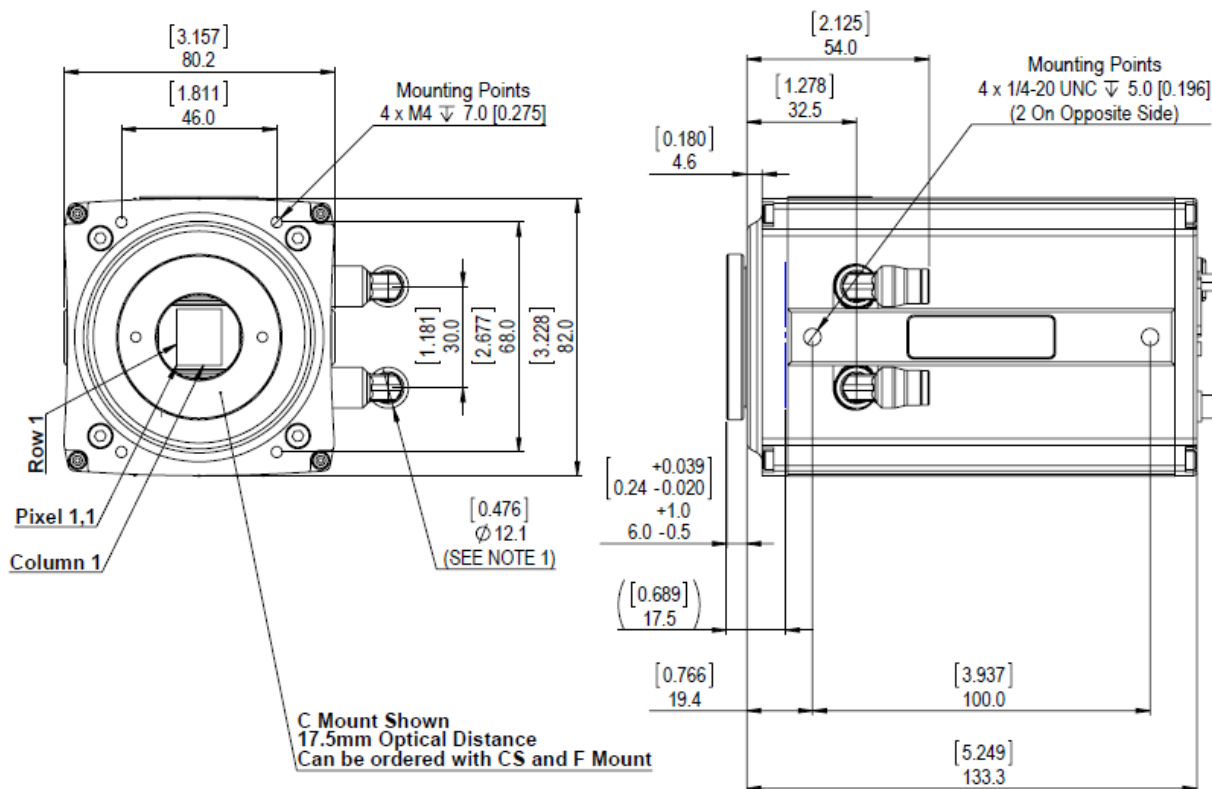
Below: Zyla Mechanical Drawings for water cooled versions of Zyla 5.5 and Zyla 4.2.



NOTES

1. WATER TAPS ACCEPT 6MM WATER PIPE. 2 X WATER ADAPTERS ALSO SUPPLIED TO ACCEPT 1/4" HOSE
2. PLEASE LEAVE 70mm MINIMUM FOR THE BEND RADIUS OF THE USB3 CABLE

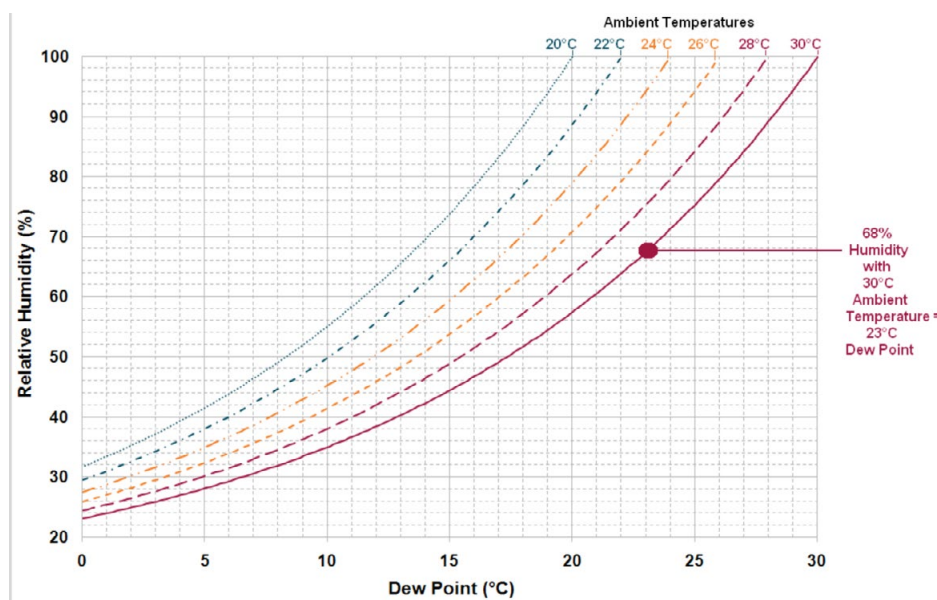APPENDIX B: DEW POINT GRAPH



The relationship between Relative Humidity and Dew Point at varying Ambient Temperature is shown below. This can be used to calculate the minimum temperature the cooling water should be set to.

APPENDIX C: OTHER INFORMATION

C.1 Terms and Conditions of Sale and Warranty Information

The terms and conditions of sale, including warranty conditions, will have been made available during the ordering process. The current version may be viewed at: http://www.andor.com/pdfs/literature/Andor_Standard_Warranty.pdf

C.2 Waste Electronic and Electrical Equipment Regulations 2006 (WEEE)

The company's statement on the disposal of WEEE can be found in the Terms and Conditions.

